



SERVOPRO Chroma Trace Gas Analyser

User Manual

Part Number: Revision: Language: 04400/001 2 UK English



This page intentionally blank

TABLE OF CONTENT

Section

Page

1	Descrip	tion and Definitions	8
	1.1 Sc	ope of this manual	8
	1.2 Sa	fety information	8
	1.3 De	scription	9
	1.3.1	Servopro Chroma Complete Rear Connector	9
	1.3.2	Servopro Chroma Reduced Rear Connector	10
	1.3.3	FID Modification	11
	1.3.4	Auxiliary Oven Rear Connecton	11
	1.3.5	Standalone PC Rear Connecton	12
2	Softwar	e license	14
	2.1 Mi	crosoft Software License Term for Windows XP Embedded	14
3	Specific	cations	19
	3.1 Ge	neral	19
	3.2 Ga	S	21
	3.3 Ov	en Specification	22
	3.4 En	vironmental Limits	22
	3.5 Di	gital Relay Outputs	22
	3.6 4-2	20 mA Outputs	23
	3.7 Di	gital Inputs	23
	3.8 Hy	drogen Relay Output	23
	3.9 Etl	nernet Port	24
	3.10 RS	-232 Port	24
	3.11 RS	-485 Port	24
	3.12 US	B Connector	24
	3.13 Ma	ains input cables (Master/Slave/Stand alone PC)	25
	3.14 Sta	indard features	25
4	System	Description	26
	4.1 Int	roduction	26
	4.2 Pla	isma cell	27
	4.3 Ele	ectric and electronic hardware description	29
	4.3.1	Signal conditioning	29
	4.3.2	Main PC and graphic display	29
	4.3.3	Oven heater and control	29
	4.3.4	I/O board	30
5	Unpack	the Servopro Chroma	31
6	Servopi	o Chroma User Interface	32
	6.1 Us	er interface control description	32
	6.1.1	Edit Box control	32
	6.1.2	Radio control	32
	6.1.3	Slide control	33
	6.1.4	Push button control	33
	6.1.5	Grid control	34

6.1.6	Combo control	34
6.1.7	Check Box control	35
6.1.8	Tab control	35
6.2 Mei	nu Description	36
6.2.1	Run	36
6.2.1.1	Real Time Chromatogram	36
6.2.	1.1.1 Historic values menu	40
6.2.	1.1.2 Alarm historic menu	41
6.2.2	Diagnostic	44
6.2.2.1	Diagnostic	44
6.2.2.2	2 Analyse Chromatogram	47
6.2.2.3	3 Trending	50
6.2.2.4	Advanced Diagnostic	53
6.2.3	Configuration	55
6.2.3.1	System	55
6.2.3.2	2 PID	59
6.2.3.3	B Flow	. 59
6234	Cycle	63
6234	5 Date and Time	69
6236	Advanced Configuration	70
6237	7 Alarm	
6238	Printers	
624	Calibration	/ / 78
6241	Calibration	70
624	Calibration of the 4.20mA outputs	/0 Q1
6 2 5	Pomoto	01
6251	Depart Software	05
0.2.3.	Demote Control	03
0.2.3.2	A heavet	84
0.2.0	About	83
6.2.7		80
6.2.8	Quit	8/
/ Installati	on and Setup	88
7.1 Sys	tem power up	88
7.2 Elec	ctrical installation	88
7.2.1	Electrical safety	88
7.2.2	Master chassis connection	90
7.2.3	Slave chassis connection	91
7.2.4	Stand alone PC with slave chassis	92
7.3 Cor	necting Servopro Chroma rear connector	93
7.3.1	Digital input	93
7.3.2	Digital output contacts	94
7.3.3	4-20 mA outputs	95
7.3.4	FID fuel shut-off relay/Hydrogen relay output	96
7.3.5	RS-485 port	97
7.3.6	RS-232	97
7.3.7	Auxiliary oven port	97

7.4 Gas installation	98
7.4.1 Introduction	99
7.4.2 Carrier Gas Hardware Selection	100
7.4.3 Connecting the Servopro Chroma	101
8 General Operation	102
8.1 Routine maintenance	102
9 Storage and Disposal	103
9.1 Storage	103
9.2 Disposal	103
10 Compliance	104
11 FID Addendum	105
11.1 Introduction	105
11.2 Overview	105
11.3 Logarithm based amplifier	105
11.4 MENUS	106
11.4.1 CONFIGURATION-ADVANCE MENU	106
11.4.2 CONFIGURATION-ADVANCE FID MENU	107
11.4.3 CONFIGURATION-ADVANCE DETECTOR MENU	109
11.4.4 CONFIGURATION-ADVANCE FLOW MENU	110
11.4.5 CONFIGURATION-ADVANCE OVEN MENU	111
11.4.6 MONITORING AND CONFIGURATION MENUS	112
11.4.7 DIAGNOSTIC-DIAGNOSTIC MENU	115
11.4.8 CONFIGURATION-CYCLE MENU	116
11.5 Startup	117
11.6 System errors	118
11.7 Troubleshooting	120
12 RS-232 Addendum	121
12.1 Start analysis	121
12.2 Analyser status	121
12.3 Send results	122
12.4 Enable-Disable Calibration	123
12.5 Respan	123
13 Product disposal In accordance with the Waste Electrical and Electronic	
Equipment (WEEE) Directive	124
14 REACH Regulation	124

LIST OF TABLES

Table name

Page

Table 1 : General Specification	. 21
Table 2 : Gas Specification	21
Table 3 : Oven specification	. 22
Table 4 : Environmental limits	. 22
Table 5 : Digital relay outputs	. 22
Table 6 : 4-20 mA outputs	. 23
Table 7 : Digital inputs	. 23
Table 8 : Hydrogen relay output	. 24
Table 9 : Ethernet cable	. 24
Table 10 : RS-232 port	. 24
Table 11 : RS-485 port	. 24
Table 12 : RS-485 port	. 24
Table 13 : Mains cable	. 25
Table 14 : Run time Menu	. 40
Table 15 : Diagnostic menu	. 46
Table 16 : Diagnostic menu 2	. 46
Table 17 : Configuration menu	. 49
Table 18 : Trending menu	. 52
Table 19 : Unregulated flow	. 57
Table 20 : Oven set point	. 57
Table 21 : Flow Factor	61
Table 22 : Peak Configuration Columns	65
Table 23 : Peak Configuration Menu	65
Table 24 : Digital Output Contacts	. 94
Table 25 : 4-20 mA connections and associated fuses	.95
Table 26 : Advanced FID menu parameters 1	107
Table 27 : Gas flow set point	117
Table 28 : Gas flow set point 2 1	117
Table 29 : Gas ignition Point 1	117
Table 30 : Troubleshooting 1	120
Table 31 : RS-232 Start Analysis	121
Table 32 : RS-232 Analyser Status 1	121
Table 33 : RS-232 Send Results 1	122
Table 34 : RS-232 Enable-Disable Calibration 1	123
Table 35 : RS-232 Respan 1	123
•	

LIST OF FIGURES

Figure name

Page

Figure 1 : Servopro chroma complete rear connection	9
Figure 2 : Servopro chroma reduced rear connection	10
Figure 3 : FID Modification	11
Figure 4 : Auxiliary oven rear connection	11
Figure 5 : Standalone PC rear connection	12
Figure 6: Edit Box control	. 32
Figure 7 : Radio control	. 32
Figure 8 : Slide control	. 33
Figure 9 : Button control	. 33
Figure 10 : Grid control	. 34
Figure 11 : Combo Box control	. 34
Figure 12 : Check Box control	35
Figure 13 : Check Box control.	35
Figure 14 : Menu description	36
Figure 15 : Real time menu	. 37
Figure 16 : Diagnostic menu	. 44
Figure 17 : Analyse chromatogram	. 47
Figure 18 : Trending menu	50
Figure 19 : Trending menu valve state	51
Figure 20 : Advanced diagnostic menu	53
Figure 21 : System menu	55
Figure 22 : PID menu	. 59
Figure 23 : Unregulated flow configuration menu	. 60
Figure 24 : Regulated flow configuration menu	61
Figure 25 : General cycle menu	63
Figure 26 : Peak configuration menu	64
Figure 27 : Peak detection configuration menu	. 66
Figure 28 : Valve configuration menu	. 67
Figure 29 : Typical sequence for valve timing	. 67
Figure 30 : Flow timing configuration menu	68
Figure 31 : Date and Time menu	69
Figure 32 : Advanced configuration menu	70
Figure 33 : Detector configuration menu	72
Figure 34 : Advanced flow configuration menu	.73
Figure 35 : Advanced valve configuration menu	.74
Figure 36 : Advanced oven configuration menu	75
Figure 37 : Alarm outputs configuration menu.	.76
Figure 38 : Printers configuration menu	77
Figure 39 : Calibration menu	78
Figure 40 : Span calibration menu	.79
Figure 41 · 4-20 mA outputs calibration menu	81
Figure 42 · Remote software menu	83
	00

Figure 43 : Remote control menu	84
Figure 44 : About menu	85
Figure 45 : Chromatogram saving menu	86
Figure 46 – Master chassis electrical connections	90
Figure 47 – Master and slave chassis electrical connections	91
Figure 48 – External PC connections	92
Figure 44– Digital input connection	93
Figure 45 – Digital outputs connection	94
Figure 46– 4-20 mA schematic and external connections	95
Figure 47 – Connecting a fuel shut-off valve	96
Figure 48: Advanced menu 1	06
Figure 49: Advanced FID menu 1	07
Figure 50: Advanced detector menu 1	09
Figure 51: Advanced flow menu	10
Figure 52: Advanced oven menu 1	11
Figure 53: Monitoring fig 1 1	12
Figure 54: Monitoring fig 2 1	12
Figure 55: Monitoring fig 3 1	13
Figure 56: Monitoring fig 4 1	13
Figure 57: Monitoring	14
Figure 58 : Diagnostic menu 1	15
Figure 59: Cycle menu	16
Figure 60 : The WEEE label 1	24
Figure 44– Digital input connection Figure 45 – Digital outputs connection Figure 46– 4-20 mA schematic and external connections Figure 47 – Connecting a fuel shut-off valve Figure 48: Advanced menu Figure 49: Advanced FID menu Figure 50: Advanced detector menu Figure 51: Advanced flow menu Figure 52: Advanced oven menu Figure 53: Monitoring fig 1 Figure 54: Monitoring fig 2 Figure 55: Monitoring fig 3 Figure 56: Monitoring fig 4 Figure 57: Monitoring fig 4 Figure 58: Diagnostic menu Figure 58: Cycle menu Figure 59: Cycle menu Figure 60: The WEEE label	93 94 95 96 106 107 109 110 112 112 113 114 115 116 124

1 DESCRIPTION AND DEFINITIONS

1.1 Scope of this manual

This manual provides installation, operation and routine maintenance instruction for the Servomex Servopro Chroma online total hydrocarbons analyser, abbreviated Servopro Chroma where applicable in the reminder of this manual.

1.2 Safety information

Read this manual and ensure that you fully understand its content before you attempt to install, use or maintain the Servopro Chroma. Important safety information is highlighted in this manual as WARNINGs and CAUTIONs, which are used as follows.



WARNING

Warnings highlight specific hazards which, if not taken into account, may result in personal injury or death.

CAUTION

Cautions highlight hazards which, if not taken into account, can result in damage to the Servopro Chroma or other equipment or property.

This manual also incorporates "Be aware of" information, which is used as follows:



This highlights information which it is useful for you to be aware of (for example, specific operating conditions, and so on).

1.3 Description



1.3.1 Servopro Chroma Complete Rear Connector

Figure 1 : Servopro chroma complete rear connection

This rear panel is present on all master units and on slave units in configurations involving a standalone pc.

- 1. Power inlet with switch and fuse
- 2. Gas trap inlet
- 3. RS-232 socket
- 4. Auxiliary oven
- 5. RS-485 socket
- 6. Gas trap outlet
- 7. Ethernet port
- 8. USB port
- 9. Analogue 4-20 mA outputs
- 10. Digital outputs
- 11. Digital input

- 12. Earth terminals for cable screens
- 13. Carrier gas inlet to detector 1
- 14. Carrier gas inlet to detector 2
- 15. Vent for carrier gasses
- 16. Vent from detector 1
- 17. Vent from detector 2
- 18. Vent from detector 3
- 19. Supporting gas inlet
- 20. Sample gas vent
- 21. Sample gas inlet to detector 1
- 22. Sample gas inlet to detector 2

Many of these ports are optional and may not be present on a particular analyser, depending on the configuration and number of detectors and intended use.



1.3.2 Servopro Chroma Reduced Rear Connector

Figure 2 : Servopro chroma reduced rear connection

This rear panel is present on all slave units in a system which contains a master unit.

- 1. Power inlet with switch and fuse
- 2. Gas trap inlet
- 3. Gas trap outlet
- 4. Auxiliary oven
- 5. Ethernet port
- 6. RS-485 port
- 7. Carrier gas inlet to detector 1
- 8. Carrier gas inlet to detector 2

- 9. Vent for carrier gasses
- 10. Vent from detector 1
- 11. Vent from detector 2
- 12. Vent from detector 3
- 13. Sample gas vent
- 14. Sample gas inlet to detector 1
- 15. Sample gas inlet to detector



1.3.3 FID Modification

Figure 3 : FID Modification

When a FID detector is installed in a Servopro Chroma, the rear connector is modified to allow for the different connections required. These modifications are:

- 1. Carrier inlet 1 is replaced by the FID vent
- 2. Detector vent 1 is replaced by a fuel inlet
- 3. Detector vent 2 is replaced by an air inlet
- 4. The supporting gas inlet is replaced by H2 contacts, which control the fuel shutoff valve (see section 7.3.5)

See section 10 for more information on the FID detector.

1.3.4 Auxiliary Oven Rear Connecton



Figure 4 : Auxiliary oven rear connection

- 1. Power inlet
- 2. Sample inlet
- 3. Sample outlet



1.3.5 Standalone PC Rear Connecton

Figure 5 : Standalone PC rear connection

- 1. Power inlet with switch and fuse
- 2. RS-232 port for slave unit 1 (marked "To Analyser")
- 3. RS-232 port for slave unit 2 (marked "Aux")
- 4. USB port
- 5. Ethernet port

CAUTION

To comply with the requirements of the EU EMC Directive, data cables used with the standalone PC (USB, Ethernet,) must be fitted with ferrite beads (Würth 74271131 or 74271221, depending on cable size), these beads should be clamped around 2 turns of cable.

CAUTION

To comply with the requirements of the EU EMC Directive, the front USB port should be use with a USB memory stick only. Rear USB port to be used with Keyboard mouse or USB memory stick only. Only keyboard supplied by Servomex shall be used

CAUTION

To comply with the requirements of the EU EMC Directive, a single turn ferrite (Wurth type 742-711-31) shall be fitted to the RS232 output cable.



WARNING

This analyser is not a medical device as defined in the medical devices directive 93/42/EEC and is not intended to be used on human beings for the diagnosis, prevention, monitoring, treatment or alleviation of disease, injury or replacement or modification of the anatomy.



WARNING

The Servopro Chroma must not be used as personal protective equipment.



WARNING

Improper installation, operation or service of this analyser may cause damage to the analyser and void the manufacturer's warranty



WARNING

All connections must be properly connected and leak free. Improper gas connection could result in explosion or death.



WARNING

Do not operate unless the cabinet is securely closed. Servicing this instrument implies possible exposure to shock hazard level voltages which cause death or serious injury.



WARNING

Any equipment connected to standalone PC shall be separated by mains voltage by at least double insulation.

The Servopro Chroma is a rack-mount type analyser to be used in fixed installations.

2 SOFTWARE LICENSE

2.1 Microsoft Software License Term for Windows XP Embedded

Microsoft Software License Terms for:

Windows XP Embedded and Windows Embedded Standard Runtime

These license terms are an agreement between you and Servomex. Please read them. They apply to the software included on this device. The software also includes any separate media on which you received the software.

The software on this device includes software licensed from Microsoft Corporation or its affiliate.

The terms also apply to any Microsoft

- Updates,
- Supplements,
- Internet-based services, and
- Support services

for this software, unless other terms accompany those items. If so, those terms apply. If you obtain updates or supplements directly from Microsoft, then Microsoft, and not Servomex, licenses those to you.

As described below, using some features also operates as your consent to the transmission of certain standard computer information for Internetbased services.

By using the software, you accept these terms. If you do not accept them, do not use or copy the software. Instead, contact *Servomex* to determine its return policy for a refund or credit.

If you comply with these license terms, you have the rights below.

1. Use Rights.

You may use the software on the device with which you acquired the software.

2. Additional Licensing Requirements and/or Use Rights.

- **a. Specific Use.** *Servomex* designed this device for a specific use. You may only use the software for that use.
- **b. Other Software.** You may use other programs with the software as long as the other programs
 - Directly support the manufacturer's specific use for the device, or
 - Provide system utilities, resource management, or anti-virus or similar protection.

Software that provides consumer or business tasks or processes may not be run on the device. This includes email, word processing, spreadsheet, database, scheduling and personal finance software. The device may use terminal services protocols to access such software running on a server.

c. Device Connections.

- You may use terminal services protocols to connect the device to another device running business task or processes software such as email, word processing, scheduling or spreadsheets.
- You may allow up to ten other devices to access the software to use
 - File Services,
 - Print Services,
 - Internet Information Services, and
 - Internet Connection Sharing and Telephony Services.

The ten connection limit applies to devices that access the software indirectly through "multiplexing" or other software or hardware that pools connections. You may use unlimited inbound connections at any time via TCP/IP.

- **3. Scope of License.** The software is licensed, not sold. This agreement only gives you some rights to use the software. *Servomex* and Microsoft reserve all other rights. Unless applicable law gives you more rights despite this limitation, you may use the software only as expressly permitted in this agreement. In doing so, you must comply with any technical limitations in the software that allow you to use it only in certain ways. For more information, see the software documentation or contact *Servomex*. Except and only to the extent permitted by applicable law despite these limitations, you may not:
 - Work around any technical limitations in the software;
 - Reverse engineer, decompile or disassemble the software;
 - Make more copies of the software than specified in this agreement;
 - Publish the software for others to copy;
 - Rent, lease or lend the software; or
 - Use the software for commercial software hosting services.

Except as expressly provided in this agreement, rights to access the software on this device do not give you any right to implement Microsoft patents or other Microsoft intellectual property in software or devices that access this device.

You may use remote access technologies in the software such as Remote Desktop to access the software remotely from another device. You are responsible for obtaining any licenses required for use of these protocols to access other software.

- Remote Boot Feature. If the *Servomex* enabled the device Remote Boot feature of the software, you may
 - (i) use the Remote Boot Installation Service (RBIS) tool only to install one copy of the software on your server and to deploy the software on licensed devices as part of the Remote Boot process; and
 - (ii) use the Remote Boot Installation Service only for deployment of the software to devices as part of the Remote Boot process; and

(iii) download the software to licensed devices and use it on them.

For more information, please refer to the device documentation or contact *Servomex*.

• Internet-Based Services. Microsoft provides Internet-based services with the software. Microsoft may change or cancel them at any time.

a. **Consent for Internet-Based Services**. The software features described below connect to Microsoft or service provider computer systems over the Internet. In some cases, you will not receive a separate notice when they connect. You may switch off these features or not use them. For more information about these features, visit

http://www.microsoft.com/windowsxp/downloads/updates/sp2/docs/priva cy.mspx.

By using these features, you consent to the transmission of this information. Microsoft does not use the information to identify or contact you.

b. **Computer Information**. The following features use Internet protocols, which send to the appropriate systems computer information, such as your Internet protocol address, the type of operating system, browser and name and version of the software you are using, and the language code of the device where you installed the software. Microsoft uses this information to make the Internet-based services available to you.

- Web Content Features. Features in the software can retrieve related content from Microsoft and provide it to you. To provide the content, these features send to Microsoft the type of operating system, name and version of the software you are using, type of browser and language code of the device where the software was installed. Examples of these features are clip art, templates, online training, online assistance and Appshelp. These features only operate when you activate them. You may choose to switch them off or not use them.
- **Digital Certificates.** The software uses digital certificates. These digital certificates confirm the identity of Internet users sending X.509 standard encrypted information. The software retrieves certificates and updates certificate revocation lists. These security features operate only when you use the Internet.
- *Auto Root Update.* The Auto Root Update feature updates the list of trusted certificate authorities. You can switch off the Auto Root Update feature.
- *Windows Media Player.* When you use Windows Media Player, it checks with Microsoft for
 - Compatible online music services in your region;
 - New versions of the player; and
 - Codecs if your device does not have the correct ones for playing content. You can switch off this feature. For more information, go to:

http://microsoft.com/windows/windowsmedia/mp10/privacy.as px.

Windows Media Digital Rights Management. Content owners use Windows Media digital rights management technology (WMDRM) to protect their intellectual property, including copyrights. This software and third party software use WMDRM to play and copy WMDRMprotected content. If the software fails to protect the content, content owners may ask Microsoft to revoke the software's ability to use WMDRM to play or copy protected content. Revocation does not affect other content. When you download licenses for protected content, you agree that Microsoft may include a revocation list with the licenses. Content owners may require you to upgrade WMDRM to access their content. Microsoft software that includes WMDRM will ask for your consent prior to the upgrade. If you decline an upgrade, you will not be able to access content that requires the upgrade. You may switch off WMDRM features that access the Internet. When these features are off, you can still play content for which you have a valid license.

c. *Misuse of Internet-based Services*. You may not use these services in any way that could harm them or impair anyone else's use of them. You may not use the services to try to gain unauthorized access to any service, data, account or network by any means.

- 4. Windows Update Agent (also known as Software Update Services). The software on the device includes Windows Update Agent ("WUA") functionality that may enable your device to connect to and access updates ("Windows Updates") from a server installed with the required server component. Without limiting any other disclaimer in this Microsoft Software License Terms or any EULA accompanying a Windows Update, you acknowledge and agree that no warranty is provided by MS, Microsoft Corporation or their affiliates with respect to any Windows Update that you install or attempt to install on your device.
- **5. Product Support.** Contact *Servomex* for support options. Refer to the support number provided with the device.
- **6. Backup Copy.** You may make one backup copy of the software. You may use it only to reinstall the software on the device.
- 7. Proof Of License. If you acquired the software on the device, or on a disc or other media, a genuine Certificate of Authenticity label with a genuine copy of the software identifies licensed software. To be valid, this label must be affixed to the device, or included on or in Servomex's software packaging. If you receive the label separately, it is not valid. You should keep the label on the device or packaging to prove that you are licensed to use the software. To identify genuine Microsoft software, see http://www.howtotell.com.
- 8. Transfer to a Third Party. You may transfer the software only with the device, the Certificate of Authenticity label, and these license terms directly to a third party. Before the transfer, that party must agree that these license terms apply to the transfer and use of the software. You may not retain any copies of the software including the backup copy.
- 9. Not Fault Tolerant. The software is not fault tolerant. Servomex installed the software on the device and is responsible for how it operates on the device.
- **10.Restricted Use.** The Microsoft software was designed for systems that do not require fail-safe performance. You may not use the Microsoft software in any device or system in which a malfunction of the software would result in

foreseeable risk of injury or death to any person. This includes operation of nuclear facilities, aircraft navigation or communication systems and air traffic control.

- 11.No Warranties for the Software. The software is provided "as is". You bear all risks of using it. Microsoft gives no express warranties, guarantees or conditions. Any warranties you receive regarding the device or the software do not originate from, and are not binding on, Microsoft or its affiliates. When allowed by your local laws, Servomex and Microsoft exclude implied warranties of merchantability, fitness for a particular purpose and non-infringement.
- 12.Liability Limitations. You can recover from Microsoft and its affiliates only direct damages up to two hundred fifty U.S. Dollars (U.S. \$250.00). You cannot recover any other damages, including consequential, lost profits, special, indirect or incidental damages.

This limitation applies to:

- Anything related to the software, services, content (including code) on third party internet sites, or third party programs; and
- Claims for breach of contract, breach of warranty, guarantee or condition, strict liability, negligence, or other tort to the extent permitted by applicable law.
- It also applies even if Microsoft should have been aware of the possibility of the damages. The above limitation may not apply to you because your country may not allow the exclusion or limitation of incidental, consequential or other damages.
- **13.Export Restrictions.** The software is subject to United States export laws and regulations. You must comply with all domestic and international export laws and regulations that apply to the software. These laws include restrictions on destinations, end users and end use. For additional information, see www.microsoft.com/exporting.

3 SPECIFICATIONS



WARNING

You must install and use the Servopro Chroma in accordance with the requirements of this section and subsequent sections of the manual. If you do not, the protection facilities incorporated into the design of the Servopro Chroma may not operate as intended, sample gas measurements may not be accurate, or the Servopro Chroma may be damaged.

CAUTION

To comply with the applicable requirements of EU EMC Directive: The case of the Auxiliary oven shall be bonded to the case of the Master/Slave unit using braid less than 20 cm long.

CAUTION

To comply with the applicable requirements of EU EMC Directive: The case of the Stand Alone PC shall be bonded to the case of the Master/Slave unit using braid less than 20 cm long.

CAUTION

To comply with the applicable requirements of EU EMC Directive: The assembly shall be bonded to a local EMC earth. The bond shall be less than 1 m in length.

CAUTION The Master, Slave, Stand alone PC and Auxiliary oven enclosures does not contain any user replaceable parts inside.

SERVOPRO CHROMA Gas Analyser

3.1 General

Master chassis dimensions: Height x width x depth:	178 x 483 x 610 mm (7 x 19 x 24 ins)
Slave chassis dimensions: Height x width x depth:	178 x 483 x 610 mm (7 x 19 x 24 ins)
Auxiliary oven dimensions: Height x width x depth:	83 x 483 x 457 mm (3.25 x 19 x 18 ins)
Stand alone PC dimensions: Height x width x depth:	178 x 483 x 130 mm (7 x 19 x 5.125 ins)
Master chassis mass:	11 - 27 kg
Slave chassis mass:	11 - 27 kg
Auxiliary oven mass:	7.5 kg
Stand alone PC mass:	5 kg
Electrical supply requirements: Voltage: Frequency:	100 to 120 Vac or 220 to 240 Vac ¹ , ² 50 to 60 Hz
Master chassis power (maximum):	400 VA
Slave chassis power (maximum):	400 VA
Auxiliary oven power (maximum):	400 VA
Stand alone PC power (maximum):	60 VA
Fuse rating/type: Master/Slave chassis Stand alone PC	Two fuses 4 amp / 250 V, Time delay, Type "T" or equivalent miniature fuses 5 x 20 mm
	Two fuses 1 amp / 250 V, Time delay miniature fuses 5 x 20 mm

 ¹ The analyser is supplied configured for operation with one of these voltage ranges. You must specify the voltage range when you order the analyser
 ² This analyser is rated in accordance to IEC 60664-3 Over Voltage Category II, Pollution Degree

Carrier gas pressure rating	
Minimum (plasma):	630.6 kPa (90 psig)
Minimum (FID):	828.6 kPa (120 psig)
Maximum:	828.6 kPa (120 psig)
Sample gas pressure rating Minimum: Maximum:	69.0 kPa (10 psig) 137.9 kPa (20 psig)
Detector type:	Plasma, TCD and FID

Table 1 : General Specification

3.2 Gas

 \sim

Carr	Type	Argon, Helium or Neon
	Pressure range	630.6 to 690.5 kPa (90 to 100 psig)
	Flow	30 ml/min to 1000 ml/min
Sam	ple gas: Type	Hydrogen, Oxygen, Nitrogen, Argon, Neon, Helium, Carbon dioxide, Carbon monoxide
	Pressure range	69.0 to 137.9 kPa (10 to 20 psig)
	Flow	25 ml/min to 150 ml/min
FID	Specific gas information ³ Gas used for combustion	Air and Hydrogen ⁴
	Pressure range for Air and Fuel	103.4 to 206.9 kPa (15 to 30 psig)
	Air flow	300 to 500 ml/min
	Fuel flow	30 to 100 ml/min
Table 2 : Gas Specification		

³ Only if FID detector is installed ⁴ Hydrocarbon traps must be used to remove hydrocarbons

3.3 Oven Specification

Chromatographic oven:		
Maximum temperature	300 °C	
Minimum temperature	Ambient	
Auxiliary oven:		
Maximum temperature	300 °C	
Minimum temperature	Ambient	
FID Oven:		
Maximum temperature	100 °C	
Minimum temperature	Ambient	
TCD oven:	200.40	
Maximum temperature	300 °C	
	A 11 /	
Minimum temperature	Ambient	
Table 3 : Oven specification		

3.4 Environmental Limits

Ambient temperature range	$+5^{\circ}C$ to $+40^{\circ}C$
Operating ambient humidity range	0 to 95% RH non-condensing
Maximum Operating altitude	2000 m (max)
Ingress protection	IP20

 Table 4 : Environmental limits

3.5 Digital Relay Outputs

Maximum voltage:	24 VDC
Maximum current:	1 Amp
Cable requirements: Type:	Multi-core cable with overall screen
Maximum size:	0.82 mm ² (18 AWG)
Maximum length	30 m

 Table 5 : Digital relay outputs

3.6 4-20 mA Outputs

3.7

Maximum load:	500 Ω
Isolation voltage:	500 V (dc or ac)
Output range Normal sample measurement Under range Over range Cable requirements: Type:	4 to 20 mA Less than 4 mA Over than 20 mA (24 mA maximum) Multi-strand twisted pair with overall screen
Maximum size:	0.82 mm ² (18 AWG)
Maximum length Table 6 : 4-20 mA outputs	30 m
Digital Inputs Maximum voltage:	24 VDC
Maximum current:	1 Amp
Cable requirements:	Multi core coble with evenall core or

Type:Multi-core cable with overall screenMaximum size:0.82 mm² (18 AWG)Maximum length30 mTable 7 : Digital inputs

3.8 Hydrogen Relay Output

This relay is only available when a FID detector is installed in the Servopro Chroma.

Maximum voltage:	24 VDC
Maximum current:	1 Amp
Cable requirements: Type:	Multi-core cable with overall screen
Maximum size:	0.82 mm ² (18 AWG)
Maximum length	30 m

Table 8 : Hydrogen relay output

3.9 Ethernet Port

Cable requirements Standard

Connector type

10 BaseT

RJ45 (CAT5 UTP)

Male 9 way D connector

Maximum length30 mCaution: This cable shall not be routed outside the building in which the analyser
is installed without the application of screening or similar protection.Table 9 : Ethernet cable

3.10RS-232 Port

Туре

Cable requirements

Maximum length Table 10 : RS-232 port

3.11 RS-485 Port

Туре

Cable requirements

Maximum length Table 11 : RS-485 port

3.12 USB Connector

Type Cable requirements

Maximum length Table 12 : RS-485 port

D-sub HD15 Male/Male shielded

Wired straight Shielded cable

NULL modem

Shielded cable

10 m

0.75 m

USB Type A Shielded standard USB cable with male Type A connector 1 m

3.13 Mains input cables (Master/Slave/Stand alone PC)

Туре	IEC 60320-1-C13
Cable requirements	-Screened
-	-Specified for 10 A / 250 Vac
	-Fit with a main plug that meets the
	requirement of the country where
	analyser us used.
Maximum length	2 m
Table 13 : Mains cable	

3.14 Standard features

- Isolated 4-20 mA output, used as process value transmission up to 8 outputs
- High resolution isolated mA chromatogram output
- Remote range identification contact output, one per peak, up to 8
- Two alarm dry contact outputs, user pre-settable limits
- One digital isolated inputs
- System status dry contact output
- Self-diagnostic system software
- Electronic carrier flow monitoring
- Electronic sample flow control system
- Colour graphic displayEthernet connection

4 SYSTEM DESCRIPTION

4.1 Introduction

The Servopro Chroma trace gas analyzer system is a basic mainframe configured for the final application. Each system is shipped pre-configured and ready to run.

The most frequent application is the measurement of impurities in bulk gases, i.e. H_2 , O_2 , N_2 , CH_4 , CO, CO_2 and NMHC in H_2 , O_2 , N_2 , Ar and He. Other applications are feasible like the sub ppb measurement of Argon

How the instrument is configured, i.e. detector, columns and valves are application dependent. The Servopro Chroma can be fitted with detector types T.C.D. (Thermal Conductivity Detector), F.I.D. (Flame Emission Detector) and Plasma Emission Detector.

When an application asks for it, a Servopro Chroma system can have one or more slave chassis to hold various detectors and columns.

A slave chassis has its own electronic system to control detector and data acquisition. All system parameters are controlled by the Servopro Chroma analytical software package.

Once properly configured, the system is started and operated as a stand alone online process analyzer with minimum operator intervention. There is no need to define any chromatographic method to detect and integrate a peak. Simple key parameters are entered into some configuration tables and it's done. The system can be calibrated and used.

Each peak is reported on screen with its process value, i.e. percentage, ppm or ppb. The software has advanced diagnostic tools and trending features that ease the trouble shooting process and configuration parameter definition.

The system comes standard with one 4-20 mA isolated output (up to 8 for the main chassis). There is also up to 8 dry contact remote range identification outputs. There are 2 operating ranges per peak with a user selectable multiplication factor between ranges of 2, 5 or 10. A high resolution isolate mA output is provided as standard feature to retransmit the real-time chromatogram signal. This output can be used with third party data acquisition on chromatographic software package. Two process alarm dry contact outputs are included. Two alarm set points per peak can be entered.

Contrôle Analytique was the first to include since the introduction of the first instrument a fail safe dry contact output used as a status alarm. This contact will be

activated in the event where there is a risk(s) that may lead to believing that the reported value may begin to be or become unreliable.

It can be seen as a common alarm for the system. All digital I/O can be configured as normally closed or open. Finally, there are 2 isolated digital inputs. One input is used as a remote start function. The other one is reserved for custom request.

The Servopro Chroma is the most complete G.C. based gas analyzer presently available.

Please, in order to appreciate and understand all the features of your Servopro Chroma system, take the time to read and understand all parts of the User's Manual.

If you have any doubt or questions as a customer of Servomex, you may contact us directly for technical support free of charge.

4.2 Plasma cell

The carrier gas flows at atmospheric pressure through a proprietary design pure quartz cell. This cell is submitted to a high frequency high intensity electromagnetic field.

The plasma cell is based on a spectroscopic emission cell, which in itself is not a new technique. On the other hand the characteristics which make this system stable and selective are the frequency, the intensity, the regulation, as well as the coupling technique and focusing (stabilizing) electrodes used to keep the plasma stable.

Under these conditions, the plasma becomes the centre of a luminous phenomenon (electroluminescence). In fact, the plasma is electromagnetically induced. Plasma is a collection of charged particles; in this case, the plasma consists of a stream of carrier gas (Ar, He or Ne). This process being an emission technique, it is very efficient for quantitative analysis. The recent advances in semiconductors as well as in optical coating and devices make the development of new instruments based on the plasma technology easier; this without the cost normally associated with this category of analytical instruments.

Once the carrier gas is ionized (charged), many spectral lines are emitted. Here, we think a few words regarding the technique used to create the plasma are necessary. There are many ways of producing light from a gas stream for analytical purposes. The electroluminescence phenomenon includes luminescence from all kinds of electrical discharges, such as sparks, arcs or tubes of different kinds, operating on direct or alternating current of low or high frequency. Some experiments were conducted in the microwave range by surface-wave induced plasma, also for the study of gas by optical emissions.

Excitation, in these cases, results mostly from electron or ion collision; that is, the kinetic energy of electrons or ions accelerated in an electric field in which the atoms or molecules of a gas are subjected to, which cause the emission of light.

By any of the above mentioned methods, characteristic emission spectra can be obtained for carrier gas and each substance in it. The emission usually varies for a given substance, depending on the mode of excitation.

4.3 Electric and electronic hardware description

4.3.1 Signal conditioning

The signal conditioning module is based on the latest technology state of the art electronic devices. Only the highest resolution with the lowest noise is used. Operational amplifiers are the best grade ones. Electrical resistors have the lowest noise and zero thermal coefficient. This leads to an analogue signal conditioning that could have a very high gain with minimum drift and noise. A special design low noise high stability analogue power supply is on board.

The signal conditioning board can accept signals from up to 8 detectors. TCD, FID and plasma emission detectors could be used. The various signals can be monitored from the diagnostic menu. The signal conditioning module has its own microcontroller to communicate with the main PC.

4.3.2 Main PC and graphic display

The main PC board runs the Servopro Chroma chromatographic software. It manages all user interface I/O and sends the information to the I/O board and the Signal Conditioning board via RS-485. Keypad and colour graphic display are directly connected to it.

4.3.3 Oven heater and control

Up to 8 ovens can be installed in the Servopro Chroma main chassis. The temperature electronic control hardware is mounted on the I/O board. There are 8 temperature control loops (PID) in the software. The temperature is measured with a RTD. The analogue to digital resolution is 24 bit (A/D converters). The heater is controlled in a pulse with modulation (PWM) scheme.

There are 8 electronic relays mounted on the I/O board. These relays turn ON at zero crossing voltage and turn OFF at zero crossing current thus eliminating EMI. Oven circuits are also protected by a fuse mounted on the I/O board.

4.3.4 I/O board

The I/O board holds all the I/O functions of the instrument. See following table.

Digital I/O	 1 isolated digital inputs 8 remote range dry contact outputs 2 alarm dry contact outputs 1 system status dry contact output 1 extra relay 10 G.C. valve contact outputs
Analogue I/O	 Soven power relays 8 oven temperature RTD input 8 isolated process 4-20 mA output 1 high resolution isolated mA output for the chromatogram

All digital inputs and outputs, analogue inputs and outputs are transient and fuse protected. The I/O board is connected to an external 40 pin I/O connector through a flat cable connector

The 40 pin I/O connector mounted on the rear panel of the instrument may be disconnected from the analyzer without the need to unscrew each wire individually

The I/O board holds the fuses for the I/O. Each fuse is socket mounted and easily replaced.

5 UNPACK THE SERVOPRO CHROMA

- 1. Remove the Servopro Chroma and any other equipment from its packing.
- 2. Remove the wrapping plastic.
- 3. Remove protective caps from inlets.
- 4. Remove protective cap from exhaust if a FID is installed.
- 5. Inspect Servopro Chroma and other items supplied, and check that they are not damaged. If any item is damaged, immediately contact Servomex or your local Servomex agent.
- 6. If you do not intend to use the Servopro Chroma immediately:
 - Refit protective caps on inlets and exhaust if a FID is installed.
 - Place the Servopro Chroma and any other equipment supplied back in its protective packing.
 - Store the Servopro Chroma as described in Section 8.

Otherwise, read section 4 (User Interface), then continue at Sections 5 onwards to install, set up, and use the Servopro Chroma.



Retain the shipping documentation and packaging for future use (for example, return of the Servopro Chroma to Servomex for servicing or repair).

CAUTION

You must remove the protective covers as specified in Steps 3 and 4 above before you use the Servopro Chroma. If you do not, you may damage the Servopro Chroma when you try to pass gases through it.

6 SERVOPRO CHROMA USER INTERFACE

All analyzer's functions are controlled through different menus of the user interface in which controls are found. The user has to interact with these controls which are explained in this section. You must become familiar with these.

6.1 User interface control description

6.1.1 Edit Box control

The Edit Box control is used to input or view a numerical value. A typical use of this control might be the input of the cycle length.

Cycle length	0	sec
Figure (+ Edi	t Day control	

Figure 6: Edit Box control

When an Edit Box control that can accept user input is active, a cursor appears in this control.

To operate an Edit Box control from the keyboard:

- > Press LEFT or RIGHT to move the cursor.
- > Press HOME to move the cursor to the beginning of text.
- > Press END to move the cursor to the end of the text.

6.1.2 Radio control

The Radio control is used to select a mode of operation. The range mode is a good example.

Range Mode
• Manual
C Automatic

Figure 7 : Radio control

To operate an Edit Box control, the mouse or the keyboard can be used. The left and right or the up and down arrows will switch between values.

6.1.3 Slide control

The slide control is used when the user may choose between items; for example, when the user has to choose between "selected peaks" or "all peaks" in the calibration menu.

All neaks	beaks	-
, imposio		

Figure 8 : Slide control

To operate a slide from the keyboard:

- > Press UP to move the slider up one position.
- > Press DOWN to move the slider down one position.
- > Press HOME to move the slider to the top of the slide.
- > Press END to move the slider to the bottom of the slide.

To operate from the mouse, move the slider to the position desired.

6.1.4 Push button control

The user clicks on a push button to trigger an action which is displayed on the button.



To operate a button from the keyboard:

➤ Use the hot key corresponding to the button.

To operate from the mouse, left click on the button.

6.1.5 Grid control

Oven Set Point
Set Point
Value
Oven
°C
°C

The grid is used to show, change or highlight data.



To operate a grid from the keyboard:

> To change a value, enter the new value and change cell.

To operate a grid with the mouse, just left-click on the desired cell to highlight it.

6.1.6 Combo control

The Combo Box control is used to choose a value between sets of predetermined values.

Acquisition time	
1 sec	•

Figure 11 : Combo Box control

To operate a Combo Box with the keyboard, press UP and Down to change between the different values.

To operate it with a mouse, left-click on the arrow and then left-click on the desired value.

6.1.7 Check Box control

The Check Box control is used to enable a feature.

🔲 Raw counts

Figure 12 : Check Box control

To operate it with the keyboard, press the space key to make it checked (enabled) or unchecked (disabled).

To operate it with the mouse, left-click in the check box to make it checked (enabled) or unchecked (disabled).

6.1.8 Tab control

The Tab control is used to switch between different pages of data. For example, a tab in the **RUN-REAL TIME CHROMATOGRAM MENU** can be selected to toggle between each conditioning boards.

All Cond. Board	Cond. Board 1
Figure 13 : Che	eck Box control

To operate it with the mouse, just click on the desired tab.
6.2 Menu Description

The analyzer's functions are regrouped in menus which are structured as in the following figure.



Figure 14 : Menu description

6.2.1 Run

6.2.1.1 Real Time Chromatogram

Pressing CTRL-R or clicking **Run** and then **Real Time Chromatogram** on the menu bar brings you to the **RUN-REAL TIME CHROMATOGRAM MENU** (fig. 10).



Figure 15 : Real time menu

The real time chromatogram menu is the main menu of the Servopro Chroma Trace Gas Analyzer software. In this menu, you can start a cycle and visualize the resulting real time chromatogram. This chromatogram takes into account all the peak data that can be configured in the **CONFIGURATION-CYCLE MENU**.

A cycle may be executed by pressing the **Start** button and stopped by using the same button that would show **Stop**. When a cycle is started, the analyser waits for the ending time of the first peak and begins its peak detection between the starting time and the ending time. If a peak is found, the peak that is detected and used for integration is shown in yellow below the raw signal coming from the detectors.

During the cycle, the progress bar rises with time. The bar starts at the left to end at the right. The cycle length can be set in the **CONFIGURATION-CYCLE MENU**. During a peak, a green circle appears at the corresponding peak of the upper grid alarm column (Al). When the peak detection process is done, the peak area is converted into the process value and displayed in the grid beside the peak name in the value field. Then, the analyzer waits for other peaks and applies the same procedure.



The cycle continues even if you exit the REAL-TIME CHROMATOGRAM MENU.

You may also choose between two injection modes using the **Injection mode** radio button. If you select **Manual**, only one cycle is executed after the Start button is pressed. If you choose **Automatic**, the analyzer restarts new cycles until you press the **Stop** button or select **Manual**.

The **Injection Status** group box let you choose if the system must inject the sample or not. When the **ON** radio button is selected, the injection will be done. If **OFF**, it will be disabled.

A radio button in the **CONFIGURATION-SYTEM MENU** let you choose between **Manual** or **Automatic** ranging. If **Manual** is selected, you may toggle between the 2 ranges by pressing the **Range** button. The range used to do the last integration of a peak or the range selected is displayed in the *R* column, next to the peak value in the grid at the left of the chromatogram.

To change a peak range, highlight the corresponding peak in the grid and press the Range button.



When Manual ranging is selected, you can't change the Range during a peak. When Automatic ranging is selected, you can never change the Range by yourself; the range selection will be done based on the result of the last peak calculations.

Note that when calibration is enabled, *Warning: Calibration enabled* will be displayed in the chromatogram display.

Another important thing about peak values is the colour displayed in the Alarm column (Al) of the peak grid. The colour is used to display alarms for each peak. The following table represents the meaning of each colour.

Colour	Meaning
White	No alarm
Yellow	Alarm 1
Red	Alarm 2

From this menu, you can also access the alarm historic menu by pressing the **Historic Alarms** button. This button will turn red when an alarm occurs, turn yellow when the menu was opened but alarms are still active, and turn green if the problem is resolved. The **ALARM HISTORIC MENU** will be explained in section 4.2.1.1.2.

A trending can be saved by pressing the **Save** button at the end of a cycle. When the save button is pressed, a dialog box appears and asks you to enter a filename. You can load this file from the **ANALYSE CHROMATOGRAM MENU**. This menu will be explained in detail later. The button is enabled only when no cycle is in progress.

The trending can also be zoomed using the mouse. Press the mouse's left button and keep it down while moving the mouse. A rectangle while appear on the screen to identify the zone to zoom. Releasing the button will apply the zoom. Pressing the **Undo Zoom** button will show the original trending without any zoom.



Remember to delete trendings that are not used anymore with the Delete button of the ANALYSE CHROMATOGRAM MENU in order to avoid filling the analyzer hard disk.



When saving a real-time chromatogram, a pop-up will advise you if the analyzer hard-drive is full. If so, delete trendings from the ANALYSE CHROMATOGRAM MENU with the delete button.

Note that a tickmark (little vertical line) is displayed on the real-time chromatogram at every starting and ending peak to help you see the integration windows.

You can view each conditioning board's trending independently by clicking on the corresponding tab at the top of the chromatogram.

Moreover, you have access to a **Print** button, which enables you to print the chromatogram. **To add a printer, see the CONFIGURATION-PRINTERS MENU description.**

A remote countdown field shows the time left when a remote starting is activated. The remote time can be set in the **CONFIGURATION-SYTEM MENU**.

If the analyzer can hold multiple configurations, the name of the current configuration will appear in the bottom right corner of this menu. Analyzer with multiple configurations is optional and this could not apply to your machine.

The following information is also displayed in the **RUN-REAL TIME CHROMATOGRAM MENU**: the carrier flow, the sample flow, the ovens temperature and the cycle time.

Push	Hot key	Function
Button		
Start/	F2	Does an injection and begins a real-time chromatogram or
Stop		stops a cycle and sets all gas chromatogram valves at their
		original position.
Range	F5	In Manual ranging only. Change the operating range.
Historic	F3	Accesses the HISTORIC VALUES MENU
Values		
TT		
Historic	F4	Accesses the ALARM HISTORIC MENU
Alarm		
Save	F6	Saves the real-time chromatogram in a file that can be
	-	opened from the ANALYSE CHROMATOGRAM
		MENU.
Print	Alt-F2	Sends the screen to the printer.
Undo	F8	Undoes the zoom on the trending.
Zoom		

 Table 14 : Run time Menu

6.2.1.1.1 Historic values menu

Displays the last 25 measures of concentration for the selected impurity. You first have to select a peak in the upper-left grid of the **REAL-TIME CHROMATOGRAM MENU** and then to press the Historic Values (F3) button to see these values.

6.2.1.1.2 Alarm historic menu

Displays the last 200 system alarms, peak value alarms or event logs that occurred and all the current active alarms. **See HISTORIC-ALARM MENU** description for more details.

System alarms:

➤ Low sample flow:

When the sample flow goes below 10 CC/min with a sample flow set point bigger than 10 CC/min;

Low carrier flow:

When the carrier flow falls below 5 CC/min for the plasma and 2 CC/min for the TCD during 30 seconds, a "Plasma #X on I/O Board #X shut down" or "TCD #X on I/O Board #X shut down" alarm will be initiated after 30 seconds to protect the system by turning off the plasma or TCD.

Plasma shut down or TCD shut down:

When a "Low carrier flow" alarm remains active for 30 seconds, turns the plasma or TCD off.

> Plasma OFF:

When the cell signal counts are lower than the starting count, it indicates that the plasma is physically off.

Starting:

When a "Plasma off" alarm is active, no "Low carrier flow" alarm is active, the starting mode is automatic (see section 8.3.1 about the **SYSTEM CONFIGURATION MENU** for the **Starting mode** definition) and no cycle is progressing (real-time chromatogram stopped or between cycles), the plasma is restarted.

> Plasma ON:

When the cell signal counts go above the starting counts after a "Plasma OFF", it indicates that the plasma is physically on.

RTD problem:

When an oven temperature falls below 10 degrees Celsius.

Carrier flow deviation:

When the carrier flow is higher or lower than the carrier flow set point by 2 CC/min, the injection is still possible and the system continues to report process values.

• Oven temperature deviation:

When an oven temperature is higher or lower than the oven temperature set point by 2 degree Celsius.

Communication problem:

When the user interface (PC) can not exchange any data with an electronic board inside the analyzer.

Peak value alarms:

(no impact on the system status alarm dry contact output)

- Alarm 1 peak #:
 When the concentration of impurity is greater than the Alarm 1 for the peak.
- Alarm 2 peak #:
 When the concentration of impurity is greater than the Alarm 2 for the peak.
- ➤ Overscale peak #:

When the concentration of impurity is greater than the actual scale of the peak.

Peripheral alarms:

Some peripherals connected with the Servopro Chroma (FID detector, HPS²...) could show their alarms / errors status into the Alarm historic menu. Please, refer to their respective user's manual for more details.

When a problem arises, a new alarm is generated and the button of the **ALARM HISTORIC MENU** turns red. If the menu is opened and an alarm is still active, it turns yellow. If another alarm occurs, it turns red again. The button will turn green if all alarms are resolved.

Every system alarm will activate (or deactivate, depending the settings in the **ALARM CONFIGURATION MENU**) the system status alarm dry contact output. Peak value alarm 1 and 2 have each a separate dry contact output that is activated (or deactivated, still depending on the **ALARM CONFIGURATION MENU**) when the alarm thresholds are reached.

Some system functions such as a real-time chromatogram are disabled when alarms are active. Rectify the problems before proceeding with your analysis.

When the problem is resolved, the same alarm message is displayed, but with an "Ok" message (except for a "Plasma shut down" that is resolved with a "Starting" and for the "Plasma OFF" that is resolved with a "Plasma ON").

Example:

- If a "Low carrier flow" occurs, the following message is displayed: (date) Low carrier flow: (hour)
- When the flow problem is rectified, the following message is displayed: (date) Low carrier flow: Ok (hour)

6.2.2 Diagnostic

6.2.2.1 Diagnostic

Pressing **Diagnostic (CTRL-P)** on the menu bar and then on **Diagnostic** brings you to the DIAGNOSTIC-DIAGNOSTIC MENU (fig. 11). The system diagnostic menu can be used for troubleshooting or only for information about the system.

This menu is divided in hardware component groups:

- Acquisition boards: read signals coming from detectors; -
- IO boards: have several functionalities related to the system as toggling relays, applying polarisation, reading mechanically regulated flows, regulating oven temperatures, etc;
- Pressure regulators: regulate flows. Acq. Board Plasma #1 Acq. Board Plasma #2 Detector signal Gain: 70 Scale Factor: 1× • Count Volt Pre-Amp: 2 Polarity: Follower --Cell 3787200 1.128674 Chromatogram 14432185 4.301126 -8 Zero (F8) IO Board #1 | IO Board #2 Pressure Regulator Ovens Flows Name Count CC/min Name ٥Ċ ٠ Flow Name Count CC/min PSI 179508 Carrier1 24 oven #1 50 Sample1 3600 75.0 5.5 Carrier2 173481 14 oven #2 378 Air 15890 150.0 24.1 oven #3 50 H2 0.8 544 0.3 Detector Generator Power ши -0 ٠ Sample1 6621 50.0 10.0 Plasma # Power Temperature (°C) Sample2 8404 50.0 12.8 65 1 In: 29.6 Out: 30.2 -

Figure 16 : Diagnostic menu

Depending on the configuration of the analyzer, different parameters are displayed under the corresponding tab of a hardware component. The following table lists the data that can be found in the **DIAGNOSTIC-DIAGNOSTIC** MENU,

Data	Meaning					
Cell	The cell counts are the detector raw signal. This data is displayed in counts, from 0 to 16777215 and also in volts, from 0 to 5V.					
Chromatogram signal	The chromatogram signal represents the signal that will be displayed on the chromatogram and used to do the peak integrations. This data is displayed in counts, from 0 to 16777215 and also in volt, from 0 to 5 V.					
Active Detector	Displays the current detector. These detectors are configured Servomex using the CONFIGURATION-ADVANCED MENU .					
Gain	Displays the gain currently applied. This gain amplifies the data to be displayed on the real-time chromatogram.					
Pre-Amp	Displays the Pre-Amp Gain applied to the signal. This gain will affect not only the Chromatogram Signal, but also the cell signal. There are 4 stages predetermined in factory that can be selected.					
Offset	Displays the voltage applied at the input of the acquisition board in order to offset the signal. This is particularly used with a FID (flame ignition detector).					
Scale Factor	Displays the factor presently used, which can be x1, x2, x5 or x10. This factor corresponds to the multiplier applied to the gain when range 1 is in use. The multiplier can be set in the <i>Range1 fact</i> column in the Peak data grid of the CONFIGURATION-CYCLE MENU . This factor also corresponds to the gain factor of the DIAGNOSTIC-TRENDING MENU .					
Polarity	Displays the current polarity of the gain. If a gain of 100 and the "inverter" (negative) value is entered, the gain is -100 . If the polarity is changed to follower (positive), the gain is $+100$. This allows negative peak to be trend positively and properly integrated.					
Flows	Displays the carrier gas flows measured on this IO Board.					
Oven temperatures	Represents the ovens temperatures in Celsius degrees.					
Detector Generator Power	Indicates the power applied to the different plasmas.					

Temperature	Shows the temperature in Celsius degrees inside the chassis and outside the chassis of a selected IO Board. If there's more than one IO Board into the analyzer, only the inside temperature is shown for the other.
Pressure Regulator	Displays the Flow's Name, the Counts, the flow in CC/min and the pressure in PSI of the different electronic pressure regulators in the analyzer.

Table 15 : Diagnostic menu

The following table lists parameters that can be manually changed to diagnose a problem or simply to observe the effects of a change.

Parameter	Meaning					
Active Detector	Let you choose which detector should be seen by the system. Only detectors for the conditioning board selected with the tab control can be selected.					
Gain	Changes the system gain.					
Pre-Amp	Toggles between Pre-Amp gain stage which can be 1, 2, 3 or 4.					
Scale Factor	Toggles between factors which can be x1, x2, x5 or x10.					
Polarity	Changes the system polarity.					
Zero (F8)	Executes a zero, which means the baseline is reset to the current counts: when the zero is executed, the chromatogram cell counts are supposed to be near 83886076.					
Detector Generator Power These values can be changed to apply a different p the plasmas (in %).						

 Table 16 : Diagnostic menu 2

6.2.2.2 Analyse Chromatogram

Pressing CTRL-L or clicking on **Diagnostic** on the menu bar and then on **Analyse Chromatogram** brings you to the **DIAGNOSTIC-ANALYSE CHROMATOGRAM MENU** (fig. 12).

101,000; 8379940,0	00						
	\bigwedge	\bigwedge	\bigwedge				
667 1334	2001 266	8 3335 40	02 4669	5336 6003	6670 7337	8004 8	3671 9338
							>
Set Position #1 (2)	Position #2 (F3)	Del	ta (#2 - #1)	Gain :	x	Load (F6)
Time (sec)	Time (s	ec)	Time (sec)		Scale Factor :	×	Delete (F7)
Count	Count		Count		Polarity :	x	Undo Zoom (F8)
Set times as star	t/end (F4)	File Name : c:/l	<4000/spance	alib.rlt	etector No (F5)	x	Print (Alt-F2)

Figure 17 : Analyse chromatogram

The analyse chromatogram menu is used when you need to analyse a chromatogram and define the right starting and ending times for the peaks. When you open this menu, the software will ask you for a file that you have previously saved (it could either be a real-time chromatogram or a trending one). If you wish to change the chromatogram, you only have to load it using the **Load (F6)** button.

All real time chromatograms are saved with the .rlt extension and .trd for the trending.

A square cursor can be used to move through each line of the graph. You can see the position of the cursor on the top-left of the screen. The X axis shows time in 1/10 sec and the Y axis indicates values in counts.

You can scroll the graph with the scroll bar on the bottom of the graph.

You can zoom in the chromatogram by using the mouse. By clicking on the left button of the mouse and holding it while moving, you can zoom in the region you want to observe. Pressing the Undo Zoom (F8) button will display the original chromatogram.

When you visualize a real time chromatogram, you cannot differentiate each conditioning board, you only get one line on the graph. But with a trending, each conditioning board read in the **DIAGNOSTIC-TRENDING MENU** can be seen with a different colour. In this case, the **Gain**, the **Scale Factor**, the **Polarity** and the **Detector No** edit box will be enabled. In these boxes, you can see the data that was active at a precise time in relation with the cursor; by moving the cursor, the data changes.

The **Detector No (F5)** button is enabled only when a trending is loaded. It will move the cursor on the graph for the corresponding detector.

You can use the **Set times as start/end (F4)** button to directly modify the peak data. The times specified in the edit control for each position will be automatically reported in the **Start** and **End** cells of the Peak data grid in the **CONFIGURATION-CYCLE MENU** for the peak selected after having pushed the **Set times as start/end (F4)** button. To change the time for each position, move the cursor to the desire position on the graph and click on the corresponding button, either **Set Position #1 (F2)** or **Set Position #2 (F3).** The difference between both positions (delta) is automatically calculated.

You may also delete a chromatogram using the **Delete (F7)** button.



As the analyzer has limited hard disk space, it is a good habit to periodically delete unused trendings. Nevertheless, when saving a real-time chromatogram or a trending from the DIAGNOSTIC MENU, a pop-up will advise you if the analyzer hard-drive is full. If so, delete trendings from the ANALYSE CHROMATOGRAM MENU with the delete button.

Push Button	Hot key	Functions			
Set Position #1	F2	Set time and counts for the position #1			
Set Position #2	F3	Set time and counts for the position #2			
Undo Zoom	F8	Resizes the trending to the normal size.			
Delete	F7	Removes a trending from the memory.			
Load	F6	Opens and displays a trending file saved in memory			
Set times as start/end	F4	Changes the starting and the ending of the peak you select with the values of time specified by the cursors positions.			
Print	Alt-F2	Prints the screen.			
Detector No	F5	Move the cursor on the graph for the corresponding detector.			

 Table 17 : Configuration menu

6.2.2.3 Trending

Pressing **CTRL-T** or clicking on the menu bar **Diagnostic** and then on **Trending** brings you to the **DIAGNOSTIC-TRENDING MENU** (fig. 13).

	Þ
Start (F2) Generator Power (F3) Re-Zero (F4)	Re-Zero Det. (F5) Polarization (F9) Time : 0 / 200000 sec
Active Detector Manual Valve D/O Boards	Print (Alt-F2) Save (Alt-F3) Undo Zoom (F8)
Det.# Acq # Gain Factor Preamp (F6) (F10)	Flows Oven
-1 1 21 ×1 1	Name CC SetPnt Follow Name CC
2 1 21 x1 1	Sample 0 0 oven #1 50.0
✓3 1 21 ×1 1	Carrier 30 30 oven #2 377.7
	Carrier 30 30 oven #3 50.0
	oven #4 49.9

Figure 18 : Trending menu

The trending menu is useful when you need to find peaks and set the proper settings for each of them, i.e. peak starting and ending times, gain, generator power and frequency, etc... These values must then be entered in the peak data table.

Before starting a trending, be sure that all parameters set in the **CONFIGURATION-SYSTEM MENU** such as sample flow, carrier flow, oven temperature, etc., have stabilized. Be sure that the gain, the detectors and the generator power are configured correctly. Nevertheless, these parameters can be changed during the trending process, except for the detectors.

To start a trending, you just have to press the **Start (F2)** button. When a trending is in progress you cannot change the detector to be trend on the graph. To stop a trending, just press the same button that will show Stop instead of Start.

You can change the generator power by pressing **Generator Power (F3)** button. A Pop-Up will ask you a value between 0 and 100% for each generator in the analyzer.

You can do a Re-Zero for the detector indicated in the edit box control beside the **Re-Zero (F4)** button. To change the detector number, just press the **Re-Zero Det. (F5)** button and it will change the detector number in the edit box. To make a Re-Zero, just click on **Re-Zero (F4)** button.

The trending can also be zoomed using the mouse. Press the mouse's left button and keep it down while moving the mouse. A rectangle while appear on the screen to identify the zone to zoom. Releasing the button will apply the zoom. Pressing the **Undo Zoom (F8)** button will show the original trending without any zoom.

You can choose if you want to follow the valve table (timings for a valve to turn ON or OFF) or not in the **CONFIGURATION-FLOW MENU**. If the check box in the **Follow?** column is checked, the valve timing will be activated for that valve. You can also manually turn on and off a valve using the combo box located in the **State** column.

۲. (Þ
Start (F2)	Generato	r Pov	ver (F3)	Re-Zero (F4)	Re-Zei	ro Det. (F	=5) Pola	rization (F	9)	Time :)00 sec
Active Detect	or Man	ual Va	alve D/O	Boards	Pri	nt (Alt-F2	2) 58	ave (Alt-F:	3)	Undo Z	coom (F8)
Valve #	State		Follow ?		Flows				1 Г	Oven	
1	Off	•	✓		Name	CC	SetPnt	Follow		Name	∘⊂
2	Off	-	v	Check all	Sample	0	0			oven #1	50.0
3	Off	-	✓		Carrier	30	30			oven #2	377.7
4	Off	-	 Image: A start of the start of	Unchek all	Carrier	30	30			oven #3	50.0
										oven #4	49.9

Figure 19 : Trending menu valve state

When the Active Detector tab of the left-bottom grid is selected, all configurations for each detector are shown. In the **Det.** # column, you can select the detector that you want to trend on the graph. A check box marked means that the detector will be trend. The **Cond** column indicates the conditioning board associated with the detector. The **Gain** column shows the actual gain for this conditioning board. If you want to change it, just change the value in the grid. The **Gain Factor** column indicates the Range Factor presently used. To change it, use the **Gain Factor (F6)** button. You have to select just one cell in the row for the desired conditioning board. The **Peak Table** column let you choose if you want to follow the peak table and the **Baseline Re-Zero** column shows if the zero will be made between each peak. To make your choice, use the same procedure as for the Gain Factor with the **Peak Table** button and the **Baseline Re-Zero** button.

The **D/O Boards** tab is used to control the digital outputs of the D/O Boards (if any) into the analyzer.

Push Button	Hot key	Functions
Start/Stop	F2	Start and Stop a trending
Generator Power	F3	Set the generator power for the detector.
Re-Zero	F4	Executes a Zero.
Re-Zero Det.	F5	Change the detector to make the zero.
Print	Alt-F2	Print the graph
Save	Alt-F3	Save the trending
Undo Zoom	F8	Restore the zoom to the original trending

Table 18 : Trending menu

6.2.2.4 Advanced Diagnostic

Clicking **Diagnostic** on the menu bar and then **Advanced Diagnostic** brings you to the **DIAGNOSTIC-ADVANCED DIAGNOSTIC MENU** (fig. 15). This menu is used to trigger different components of the analyzer.



Figure 20 : Advanced diagnostic menu

1) 4-20 mA Diagnostic:

This section is used to manually control the 4 to 20 mA analogue outputs.

First select the board to diagnose by using the combo box just under the title "4-20mA Diagnostic". Enter a value in the grid from 4 to 20 for the corresponding 4-20 mA # and press **ENTER**. The analogue output will be set to this value and will stay at this value until it is changed from this menu or by normal operations from outside this menu.

2) Relay Diagnostic:

This section is used to verify if the Range relays on the I/O Boards are working properly. First select the I/O Board to diagnose by using the combo box just under the title "Relay Diagnostic". By using the combo box in the state column of the grid, you can switch the relay ON or OFF depending on the configuration in **CONFIGURATION-ADVANCED CONFIGURATION MENU**. The value is applied after an **ENTER** on the keypad or when a mouse click is performed outside the combo box.

3) Valves Diagnostic:

This section is used to verify if the valves on the I/O Boards are working properly. First select the I/O Board to diagnose by using the combo box just under the title "Valves Diagnostic". By using the combo box in the state column of the grid, you can switch the valve ON or OFF. The value is applied after an **ENTER** on the keypad or when a mouse click is performed outside the combo box.

4) Status, Alarm 1 and Alarm 2

These group boxes let you change the state of the corresponding relays of Alarm and Status for a short period of time. If the analyzer processes a new alarm or error, these relays will return to their normal state after a short moment even if you have triggered a different state. By default, these relays are on the first IO board.

6.2.3 Configuration

6.2.3.1 System

Pressing **CTRL-S** or by clicking **Configuration** on the menu bar and then **System** brings you to the **CONFIGURATION-SYSTEM MENU** (fig. 16). The system configuration menu is another very important menu. Many parameters are set there. These parameters are listed in the following paragraphs.

Analog © Tr O Ho	i output	-mA Failure Off C Low C High	Mode —	Generator S C Manual C Automa	itarting Mode - itic	Range C M	Mode anual utomatic	Auto-Start on Boot-Up Ves No	
-Flow Se	et Point Flow Name	Set Point CC/min	Value CC/min		-Oven Se	et Point	Set Point °C	Value °C	
	Sample	30.0	30.0			oven #1	50.0	50.0	
	Carrier 1-52	30.0	30.0			oven #2	30.0	30.0	
	Carrier 2-52	30.0	30.0			oven #3	50.0	50.0	
						oven #4	50.0	50.0	
Reload I	mote Starting sec Manufacturing	Enable Setting	Foreen S	iaver Enable sec	-Range C Peak & H2 O2	Configuratio Activ # on Ra Opened Opened	n ve State ange #2 d 💌	Lock Range (F2)	
Cl Cha	hange Passwoi inge Configura	tion	ctive Config	: None				Set all closed	

Figure 21 : System menu

1) Analogue output:

You can choose between two modes. **Hold** mode maintains the 4-20 mA outputs even if a new cycle is started. In **Track** mode, analogue outputs always track input gas values.

2) mA Failure mode:

In the case of a system status alarm, this control sets <u>all</u> the 4-20 mA outputs below 4mA if **LOW** is selected and higher than 20 mA if **HIGH** is selected. Selecting **OFF** will disable this function and the 4-20 mA outputs reflect signal values as in normal operation.

3) Generator Starting mode:

• **Manual:** the automatic re-start feature is disabled and you may enter any plasma power value between 0 and 100%. It is useful when troubleshooting or re-configuring the system.

• Automatic: when the cell raw counts are lower than the cell Plasma starting count value, a higher power will be applied to the cell to restart the plasma. <u>NOTE</u>: The system checks if the plasma is OFF between cycles. This feature is used only when the analyzer has a plasma cell.

4) Range Mode:

Automatic: automatically changes the range according to the result of the previous peak integration. If the result of peak integration gives a value higher than 99% of the range presently used, the analyser switches automatically on the next higher range. The next integration gives the proper result. If the result of a peak is lower than 90% of the lower range presently used, the analyser switches to this lower range. **Manual**: you may change range manually.



This process is independent for all peaks, which means that in Automatic ranging, the second peak range can stay the same even if the first peak changes range.

5) Auto-Start on Boot-Up:

When set to **Yes**, will try to start a cycle (or several cycles if The **Injection mode** has been set to **Automatic**) on the analyzer boot-up after all the alarms are resolved.

6) Flow Set Point:

Indicates the flows and the set point for each of them. There are two kinds of flow, regulated flow and unregulated flow. A regulated flow is controlled by an autonomous electronic board called Pressure Regulator Board and an unregulated flow is not controlled at all, but only read by a sensor managed by the I/O Board.

Regulated flow:

This value is used by the flow control loop. In the third column of the Flow Set Point grid, the real flow is displayed. There can be a delay between the time you specify a new set point and the time the real flow is displayed and stabilized to the specified set point. To change the set point, change the value in the second column for the corresponding flow. The range of the values that can be set corresponds to the range of the associated flow sensor table (see **CONFIGURATION-FLOW MENU**).

Unregulated flow:

This set point value is only used as a reference for flow deviation alarm. The real flow set point is determined by the pressure regulator on the back of the Servopro Chroma. If the actual flow is 2 CC/min lower or 2 CC/min higher than the reference, the flow deviation alarm is added in the ALARM HISTORIC and the contact status is opened (or closed, depending the configuration in the **CONFIGURATION-ALARM MENU**). This is, in fact, an alarm set point.

Minimum	0
Maximum	200
Resolution	1

 Table 19 : Unregulated flow

7) Oven set point:

Defines the oven temperature in Celsius degrees. This is used by the oven temperature control loop. The real temperature of the oven is displayed in the third column of the Oven Set Point grid. To change the set point, change the value in the second column for the corresponding oven. Be aware that the desired set point can take several minutes to be reached and stabilized.

Minimum	0
Maximum	300
Resolution	1

Table 20 : Oven set point

8) Remote Starting Enable:

When this check box is checked, a cycle can be started by the digital input #1. The edit box specifies the time that elapsed before the cycle starts after the digital input #1 is activated. When the 24 Volts is applied to the digital input #1, the remote starting countdown starts (it can be viewed in the **RUN-REAL TIME CHROMATOGRAM MENU**). To start a cycle, the voltage must not be applied anymore when the countdown reaches 0. If voltage still applied, the countdown restarts.

9) Screen Saver Enable:

When this check box is checked, the screen saver is enabled. The edit box specified the time that elapsed before the screen power off. By moving the mouse or pressing a key, the screen will power on.

10) Reload manufacturer settings:

To reload the analyzer original configuration that was done in factory, click on this button. Before clicking on this button, please, call Servomex for more information.

<u>11) Range Configuration:</u>

This is used to configure the relays on the I/O Board for the Peak Ranges. By changing the value of the combo box for the corresponding peak, you can change its state when range 2 is active.

By the Lock Range button, you can also Lock or Unlock the analyzer ranges. If you press the lock button, the range will be locked until you press the unlock button (the lock button and the unlock button are the same. Only the text changes). *The range can be lock if the analyzer is in autoranging mode, but it will be ignored*. By locking the ranges, you disable the range button in the RUN-REAL TIME CHROMATOGRAM MENU so that the ranges cannot be changed even if manual ranging is ON.

12) Print configuration

This button is used to print the system configuration. In a multiple configurations environment, this will only print the current system configuration

13) Change password

This button is used to set a new password for the Servopro Chroma. This password is used when you want to quit the software or access the **CONFIGURATION-ADVANCED MENU**.

14) Change configuration

If you buy a Servopro Chroma with more then one configuration, you can change the active configuration by clicking the "Change configuration" button. When you change the active configuration, all parameters are uploaded according to the new selected configuration.

15) Change password

By default, the general password used in the analyzer is its serial number. If, for any reason, you need to change it, click on this button. A dialog window will ask you for the old password and will require you to input the new password twice.

6.2.3.2 PID

Clicking on **Configuration** on the menu bar and then on **PID** brings you to the **CONFIGURATION-PID MENU**.

Flow	Oven
Sample 💌	1 💌
Proportional 0.02	Proportional 7.00
Integrator 0.03	Integrator 1.00
Differential	Differential 0.00
Save	Save

Figure 22 : PID menu

This menu lets you set the PID settings for flows and ovens. *Be sure to save your settings before changing Oven # and Flow #*.

6.2.3.3 Flow

Clicking on **Configuration** on the menu bar and then on **Flow** brings you to the **CONFIGURATION-FLOW MENU** (fig. 18).

This menu contains the flow sensor table. This data is set up by Servomex. If you have to recalibrate some of these data, follow these steps:

Unregulated flow:

First of all, you need to choose the flow that you want to calibrate with the combo box on the top-left of the view. There are two kind of gas: Regulated flow and unregulated flow. The Simple flow is changed using a gas regulator. For calibrating these sensors, you need a 0 CC/min point, a point higher than 200 CC/min and eight points between these values. Now, turn the gas valve and

note the Flow meter counts in the **Flow Setting Tool**. Use a bubble flow meter to find the reading of the flow in CC and write down these two values on paper. When you have 10 reading points, you must enter these in the flow table using the keyboard.

		1
Flow	Counts	-
0.0	124609	
6.4	134443	
13.6	148618	Flow Setting tools
28.2	168453	
32.9	175482	
38.2	182769	
44.7	192787	
50.3	198273	
57.9	207309	
72.6	221985	
		Flow meter counts : 0
		Flow Factor : 1.0
Validate (Fi	3) Print (Alt-F3)	1

Figure 23 : Unregulated flow configuration menu

Regulated flow:

For the regulated flows, the same principle as in the unregulated flow sensor calibration is used, but you can control the gas pressure with the controlled valve using the **Flow Setting Tool**. Wait for the PSI of the pressure valve to stabilize to the new set point. Now, use a bubble flow meter to find the reading of the flow in CC. When you have 16 reading points, you must enter these in the flow table using the keyboard. Note that the Flow Control Type depends on the tubing configuration of the system (bypass or direct). Therefore, keep the Flow Control Type as it is.

Flow	PSI	7	
0.0	1.0		
28.0	1.5		
40.0	2.0	Flow Setting tools	_
52.6	2.5		
64.9	3.0	7.0 PSI	
75.5	3.5		Flow Control Type
85.1	4.0		C Direct
93.8	4.5		• Direct
100.8	5.0		Bypass
110.1	5.5		
123.7	6.0	PSI: 1.0	
131.9	6.6		
131.9	6.6		
131.9	6.6	Actual control output: 0.0 %	
131.9	6.6	Flow value (CC) : 0.0	
131.9	6.6		
		Factor : 1.0 Slope : 986835190	

Figure 24 : Regulated flow configuration menu

In this menu, the control loops (flow set point values) are disabled. But as soon as this menu is closed, the control loops are re-enabled.

A flow association must be configured in the **CONFIGURATION-ADVANCED MENU** to know which flow is a regulated flow and which is a unregulated flow.

Flow Factor:

Many different types of gases could be handled by the Servopro Chroma. Flow transducers use the thermal conductivity principle to determine the volumetric flow (CC/min) of the gas. Various gases have different conductivities, for example, Helium and Hydrogen have a much higher thermal conductivity compared to Argon or Oxygen. The **Flow Factor** can be applied to the flow table in order to obtain the proper flow for various gases. Here are some factors for various gases:

Gas	Factor
Argon	1.0
Hydrogen	8.7
Helium	8.5
Oxygen	0.92
Nitrogen	1.04

Table 21 : Flow Factor



The minimum pressure required for proper operation could need to be re-adjusted in order to avoid oscillation of the flow control loop when changing the sample gas type. This is application dependent.



The sample flow factor can be determined by using a bubble flow meter on the sample vent. To find a new flow factor for new sample gas, you m ay use a bubble flow meter connected on the sample vent bulkhead on the rear panel of the analyzer. Enter an arbitrary factor until the sample flow displayed on the LCD is close or equal to the flow measured with the bubble flow meter, that's all! For the sample flow, the absolute value is not critical, the stability is. This means that at the end of a cycle, the sample flow must be the same as the previous one. This is important for stability.

6.2.3.4 Cycle

Clicking on **Configuration** on the menu bar and then on **Cycle** brings you to the **CONFIGURATION-CYCLE MENU** (fig. 20).

This menu contains all the information that the system needs when a cycle (Real time chromatogram) is in progress.

1) General (Cycle Data and 4-20 mA Peak Association):

This menu is used to set the duration of the cycle (length of one chromatogram) and the duration displayed in the **REAL-TIME CHROMATOGRAM MENU**. Also, it is used to assign a peak to a particular 4-20 mA output.



Don't change these values unless you know exactly what you are doing!

General			ß		
Peak	Cycle Data				
Configuration		Cycle len	gth: 500	sec	
Detection	Second o	n chromato	gram: 1000	sec	
Valve					
Flow	-4-20 mA Peak Association	on			
		Peak #	I/O Board #	4-20 mA #	
		1	1	1	▼
		2	1	1	
		3	1	1	
		4	1	1	
		5	1	1	<u>-</u>
	Save Beak As	cociation Cr	id (52)		Print (E2)
		sociación di			

Figure 25 : General cycle menu

2) Peak:

Peak Configuration:

The data that is found in the next menu is used during an analysis. The system parameters change in real time according to each peak configuration.

	Name	Avr	Start	End	Det		Alarm 1	Alarm 2	Range 2	Range 1	Un	it
Peak	Ar	1	237.0	276.0	3	•	1000000.0	1000000.0	1000	1×	▼ PPI	3 💌
Configuratio	02	1	303.0	423.0	2	•	1000000.0	1000000.0	1000	1x	💌 PPI	3
Configuration	N2	1	435.0	482.0	8	•	1000000.0	1000000.0	1000	1×	💌 PPI	M 💽
Detection	C02	1	595.0	662.0	6	•	1000000.0	1000000.0	1000	1x	💌 PPI	1
/alves	co	1	675.0	789.0	6	•	1000000.0	1000000.0	1000	1×	▼ %	•
low												
icripts	-Peak	#1 da	ita Gain: [30			Plasma	a power:	65	%		
		Po	olarity:	Follower	•		Pre-Amp	olification:	1	•		

Figure 26 : Peak configuration menu

In the grid at the top of this menu, you found general information that does not depend upon the kind of detector. Here's a description.

Column	Explanation
Name	Name of the peak.
Avr	Average that can be used to calculate the impurity value of the peak. An
	average of 1 disables the feature.
Start	The starting time of a peak.
End	The Ending time of a peak.
Det.	The detector number used to trend the peak. Refer to the CONFIGURATION-ADVANCED MENU for the detector number association.
Alarm 1	When the peak exceeds this value, the alarm 1 will be activated.

Alarm 2	When the peak exceeds this value, the alarm 2 will be activated.
Range 2	(Range 2 Scale) This value indicates the full scale of the analyzer
	(Range 2).
Range 1	(Range 1 factor) The multiplier applied to the system, when the analyzer
	is on the first range.
Unit	Unit of the peak (PPM, PPB or %)

Table 22 : Peak Configuration Columns

Underneath the grid, you find extra information about the peak that is currently selected in the grid. This information depends on the type of detector.

Polarity	The polarity: Follower or Inverter.
Gain	The peak gain.
Plasma/generator Pwr	Detector Generator Power used for the corresponding peak.
Pre-Amp	Shows the stage of the Pre-Amp for the corresponding peak.

Push Button	Hot key	Functions
Add Peak	F2	Add a peak to the grid
Delete selected peak	F3	Delete the peak selected in the grid.
Save peak table	F4	Save the peak grid
Print Peak Table	Alt-F2	Print the peak grid
Peak Detection Cfg.	None	Brings the Peak Detection Configuration Menu
T.L. 32 . D L (7 P	N/

 Table 23 : Peak Configuration Menu

Peak Detection:

In this menu, you can modify constant that are used to detect if and when there is a peak.

General						
Peak						
Configuration						
Detection						
Valve	Peak	H2	02	N2	CH4	со
	RMS Segme	4	4	4	4	4
Flow	PD segment	5	5	5	5	5
	PD Nb Segm	5	5	5	5	5
	Cut-Off fre	0.10	0.10	0.10	0.10	0.10
	Filter Length	40	40	40	40	40
	THRSH Factor	1	1	1	1	1
	LDL	0.00	0.00	0.00	0.00	0.00
	1 Derivative	1500	1500	1500	1500	1500
	2 Derivative	300	300	300	300	300
			Sav	e data		

Figure 27 : Peak detection configuration menu



AVOID CHANGING THESE SETTINGS: they are configured in factory by Servomex.

3) Valve data:

General						
Peak	Valv	e timing configuration	on			1
Configuration		1			-	
Detection		,				
Valve		0	n (sec)	Off (sec)		
Flow		0		30		
FIUW		7.	199	7200		
		7.	199	7200		
		7:	199	7200		
		7:	199	7200		
		7:	199	7200		
		7:	199	7200		
		7:	199	7200		
		7:	199	7200		
		73	199	7200		
		Save Valve #1	l timing	Р	rint (F3)	
	R					

Figure 28 : Valve configuration menu

This group box is used to control the valve state. By clicking in the combo box, you can change the time at which the valve go ON (open) or OFF (close). Each row in the grid indicates a sequence for the selected valve.



Don't forget to save valve timing data with the **Save Valve** button each time you select a new valve in the combo or if you quit the menu.



A valve timing won't be applied unless the timings are saved first.

<u>4) Flow</u>

This group box is used to configure flow timing table. A flow timing table can be used to change a flow set point during a cycle.

General		
Peak Flow timing confi	guration	
Configuration	- Carrier #2 PB #	3
Detection		
/alve	Time (sec)	CC/min
	0	30.000000
)W	7200	0.000000
	7200	0.000000
	7200	0.000000
	7200	0.000000
	7200	0.000000
	7200	0.000000
	7200	0.000000
	7200	0.000000
	7200	0.000000
	Save flo	ow timing (F2)

Figure 30 : Flow timing configuration menu

To set a flow timing table, select the proper flow name using the combo box. Then enter times and the flows into the table. A time of 7200 sec means that this point isn't use. According to figure 25, the set point will be set to 30 cc/min at the beginning of the cycle and won't change during the cycle because other timings are set to 7200 sec. Once you have fill up the grid, push the **Save flow timing** button to save data.

6.2.3.5 Date and Time

Clicking on **Configuration** on the menu bar and then on **Date and Time** brings you to the **CONFIGURATION-DATE AND TIME MENU** (fig. 26).

Date and Time	•	16:53:11	
2007-08-06	T	<u></u> 2007-08-06	
	Apply		

Figure 31 : Date and Time menu

This menu lets you set the date and time of your analyzer. By using the combo box control, you can change the value of the date and of the time. Just click on Apply to set the desired value.

6.2.3.6 Advanced Configuration

Clicking on **Configuration** on the menu bar and then on **Advanced** brings you to the **CONFIGURATION-ADVANCED MENU** (fig. 27).

This menu is the analyzer core. It defines all the hardware components on which the Servopro Chroma is built.



Therefore, you normally don't have to change these configurations; they are set in factory and reflect the hardware of the analyzer.

Before entering the Advanced Configuration menu, the analyzer will ask you for the general password of this analyzer. This password is, by default, the serial number of the analyzer, but can be changed into the CONFIGURATION-SYSTEM CONFIGURATION MENU.

If for any reason you have to change any parameters of this menu, please contact Servomex.

General Detector	General				
1					
2	Item	Quantity	System Alarm		
3	Cond. Board	1	● On ● Off		
Flow	FID	0			
FIUW	I/O Board	1	RS-232 Communication —		
Valve	Flow	3	O On 💿 Off		
Oven	Valve	4			
	Detector	3	Take Snapshot		
	Oven	4	· · · · · · · · · · · · · · · · · · ·		
	Plasma	1			
	Press. Board	3			
	Temp. Ctrl.	0			
	D/O Board	0			

1) General:



In this menu, all Servopro Chroma alarm and error monitoring can be turned off when a user wants to work with the analyzer even if for example a flow is too low for a proper operation.

<u>NOTE</u>: Turning off System Alarm can be dangerous, for example if you are working with Hydrogen: the Servopro Chroma will stop advising you of any further problems!

Also, this is where the number of electronic boards, components and functionalities are entered.

Moreover, the RS232 communication can be turned ON and OFF if the system is purchased with this option.



AVOID CHANGING THESE SETTINGS: they are configured in factory by Servomex.

Finally, the **Take Snapshot** button can be used to help Servomex diagnose a problem into your analyzer. Use this button only under recommendation from Servomex.
General	
Detector	Detector #1
1 2	
3	
5	Detector Type
6	Plasma
7	Cond Board # Channel # Location Placma #
	1 ▼ 1 ▼ 1 ▼ 1
9	
Flow	
Valve	
Oven	Save Cancel

2) Detector Association:

Figure 33 : Detector configuration menu

When using a plasma cell, a conditioning board must also be selected with the right channel. Moreover, a plasma detector must be associated to the cell by selecting its location on an IO board and its number on this IO board. The plasma detector association is used to process alarms such as Plasma Off, Plasma On and Plasma Starting.

This list includes all detectors, including those contained within any slave units connected to the master unit or standalone PC. Bear in mind however that each individual conditioning board can only process up to 8 detectors.

Don't forget to save the new configuration by clicking on the Save button.



AVOID CHANGING THESE SETTINGS: they are configured in factory by Servomex.



General Detector			Flo	w			
1							
2							
3	Name	Location		Sensor #		Associated Pla	isma
	Carrier1	I/O Board #1	•	Sensor #4	-	Plasma #1	-
5	Carrier2	I/O Board #1	-	Sensor #5	•	Plasma #2	-
6	Sample1	Press. Reg. #1	•				-
	Air	Press. Reg. #2	-				-
(H2	Press. Reg. #3	-				-
8	Sample1	Press. Reg. #4	-				-
9	Sample2	Press. Reg. #5	-				-
Flow	Carrier3	I/O Board #2	-	Sensor #3	•	Plasma #1	-
1 IOW	Carrier4	I/O Board #2	•	Sensor #4	-	FID	-
Valve							
Oven		Save		Cance	<u> </u>		

Figure 34 : Advanced flow configuration menu

The Flow menu associates a flow with a sensor on a certain I/O Board or Pressure Board. Name the flow, choose its location on the appropriate I/O Board, select Sensor number 3, 4 or 5 for a carrier, and associate a plasma detector (used to generate alarms and errors). If the location of the flow is Pressure Board, no sensor or plasma association is needed.

Don't forget to save the new configuration by clicking on the Save button.



AVOID CHANGING THESE SETTINGS: they are configured in factory by Servomex.



4) Valve:

Figure 35 : Advanced valve configuration menu

The Valve menu associates a valve number to a valve channel on an I/O Board.

General Detector	Oven
1 2	
3	Name Location Channel # Max ºC Min ºC
	Oven1 I/O Board #1 🔽 Channel #1 💌 300.00 0.00
5	Oven2 I/O Board #1 🔽 Channel #2 💌 300.00 0.00
	Oven3 I/O Board #2 🔽 Channel #1 💌 300.00 0.00
b	Oven4 I/O Board #2 🔽 Channel #2 💌 300.00 0.00
7	FID FID #1 Channel #1 300.00 0.00
8	
9	
Flow	
Val∨e	
Oven	SaveCancel

5) Oven:

Figure 36 : Advanced oven configuration menu

The Oven menu associates an oven with a certain I/O Board and a certain channel on this IO board. It is also possible to enter the maximum and minimum value for the temperature set point of this oven.

Don't forget to save the new configuration by clicking on the Save button.



AVOID CHANGING THESE SETTINGS: they are configured in factory by Servomex.

6.2.3.7 Alarm

Clicking on **Configuration** on the menu bar and then on **Alarm** brings you to the **CONFIGURATION-ALARM MENU** (fig. 32).

Status Alarn	1		
C Opena	d when acti	ve 💽 Closed when a	tive
Alarma Chata			
-Alarm State			
I/O Bo	ard #	Active State on alarm 1	Active State on alarm 2
1		Opened 🗾	Opened 🗾
	Set	All Opened	Set All Closed
	_		

Figure 37 : Alarm outputs configuration menu

This menu lets you configure the alarms. The status alarm relay can be closed or opened when a status alarm is activated. Click on the desired radio box button.

To set the active state for the Alarm 1 and Alarm 2 relay on the I/O Board, use the corresponding combo box in the Alarm State grid.

6.2.3.8 Printers

Clicking on Configuration on the menu bar and then on Printers brings you t	0
the PRINTER CONFIGURATION MENU (fig. 33).	

\\ATLAS\HP Color LaserJet 3500 \\ATLAS\HP LaserJet 5100 Series*
Add Printer Delete Printer Set as Default Printer Printer's Properties
(* = Default Printer)

Figure 38 : Printers configuration menu

The printers already installed on your analyzer will be listed in this menu. It is possible to configure a new printer by adding it with the **Add Printer** button. This will show a new dialog wizard window that will help you to choose a new printer shared over your network. The **Set as default** button determines which printer is used as default for printing. The printer's properties can be modified with the **Printer's Properties** button. To remove a printer, use the **Delete Printer** button.

6.2.4 Calibration

6.2.4.1 Calibration

By clicking **Calibration** on the menu bar and then **Calibration**, you access the **CALIBRATION-CALIBRATION MENU** (fig. 34).

Name	THRSH	Span Ga	as	Result	Rng	Select		TRS Done	Span Done	Response Factor	Area Count
N2	1352.7	21.90	-	38.63		NO	•	YES	YES	0	10644726
CO2	2224.8	102.00	-	95.72		YES	•	YES	YES	0	54307524
AR	3329.6	16.00	-	< 1.00		NO	•	YES	YES	0	2135238
02	4987.8	16.50	-	2.69		NO	•	YES	YES	0	9537891
						Tin	ne:	0 sec			
					Enable	Tin e (F2)	ne :	0 sec	Add	N2 calib. values	Enable (F3)
-Noise	Threshold Ca	alibration			Enable	Tin e (F2)	ne :	0 sec Span Calib	ration	N2 calib, values	Enable (F3)
-Noise	Threshold Ca	alibration :cted pea	ks		Enable	Tin ≥ (F2)	ne :	0 sec Span Calib 	ration Add	N2 calib. values	Enable (F3) Automatic Re-Span
-Noise	Threshold Co Sele All p	alibration :cted peal ;eaks	ks		Enable	Tin e (F2)	ne :	0 sec Span Calib 	ration Add Selected (All peaks	N2 calib. values	Enable (F3) Automatic Re-Span
-Noise	Threshold Ca Sele All p	alibration acted peal aeaks	ks		Enable	Tin e (F2)	ne :	0 sec Span Calib	ration Add Selected (All peaks	N2 calib. values	Enable (F3) Automatic Re-Span
-Noise T	Threshold Co Sele All p Start (F	alibration icted pea ieaks 4)	ks	Calculate TI	Enable	Tin 2 (F2)	ne :	0 sec Span Calib 	ration Add Selected (All peaks Start (F6)	N2 calib. values	Enable (F3) Automatic Re-Span

Figure 39 : Calibration menu

This menu is used to calibrate the system for the different gases you have to analyze.

Be sure to first configure all parameters of the system that you find in the **CONFIGURATION-SYSTEM MENU** and in the **CONFIGURATION-CYCLE MENU**. The analyzer must be properly started and purged before the execution of a calibration.

1) Noise Threshold (THRSH) Calibration:

This calibration step will find the noise level used for the peak detection process. In this step, **no injection** will be done. It is just a noise baseline on range #1, as the highest gain applied is always on this range. When starting the cycle, it will automatically switch to range #1. The RMS (Root Mean Square) value of this noise will be saved as noise and used for the peak pre-recognition. First, you need to enable the Noise Threshold Calibration with the corresponding **Enable** button. After a few cycles in automatic injection, you can press **Calculate Threshold** button. The RMS value will be shown in the THRSH column of the grid. You can **Calculate Threshold** for **All peaks** or only the **Selected peaks** by changing the corresponding slider state. When **Selected peaks** is selected, only peaks that have a **YES** in the **Select** column of the grid will be considered.

When the Noise Threshold Calibration is done, **YES** is displayed in the **TRS done** column of the grid.

2) Span calibration:

Calibration gas values must be set for each individual peak before starting a calibration run. To set calibration values, select a peak name with the mouse and then push **Add "peak name" calib values** button. A popup menu appears (Figure 35).

				,	-			1	
Name	THRSH	Span Gas	Result	Rng	Select	TRS Done	Span Done	Response Factor	Area Count
H2	1030.8	12 50 💌	20.20		NO	▼ VES	YES	0	40646428
02	356.2	Multi-point	s calibratio	n: 02		×	/ES	0	61203184
N2	48.9	Number of	f points: 2		_		/ES	0	83346720
CH4	614.2						/ES	0	111344440
co	5111.7	Index	Value	•	Are	a	/ES	0	68712560
		1	0.00		0				
		2	14.00)	542	85704			
								ĸ	
								~~~~~~~~~~~~~~~~~~~~~~~~~~~~~~~~~~~~~~~	
									(
	<b>-</b>						Add	02 calib, values	Enable (F3)
[Noise	Inreshold						ation		
	📜 Se		_	_	_		Selected	peaks 🗌	Automatic Re-Spap
						ОК			Ko-opan
	I All						All peaks		
	Start (	F4)	Calculate Th	hresho	ld (F5)		Start (F6)	RE-SPA	N (F7)

Figure 40 : Span calibration menu

First of all, you must enter the number of point that will be used to calibrate the system. Normally, user uses only two points: point 0.00 by default and another value. If you have a calibration system like Kontrol Analytik's GCS-100, you can calibrate your system for an unlimited number of impurities. Consequently,

the analyzer results will be more accurate. Once calibration values have been entered, click **OK** to go back to **Calibration** main menu. If you have more than one peak, do the same steps for the remaining peaks. Once all values are set, a calibration cycle can be performed. Click the **Enable** span calibration. Then select the impurity value for each peak using the combo box in the **span gas** column. Then, if you wish to calibrate only some peaks, select **YES** in the select column for the peaks you want to calibrate and then, select **Selected peaks** in span calibration menu. If you want to calibrate all peaks, select **All peaks**. To run a calibration cycle, push the start button and wait for cycle end. Actual cycle time is displayed under the grid. When the cycle is completed, push the **re-span** button to accept the calibration. If you want to calibrate other impurity values, select them in **span gas** column and repeat the previous steps. To disable calibration mode, click on **disable** span calibration.



*Calibration enabled* will be displayed in the **RUN-REAL TIME CHROMATOGRAM MENU**. It is suggested and of common sense to do the calibration in **Automatic** injection mode. This allows a better gas equilibrium.



When calibration is enabled, you can't change range configuration



Don't forget that prior to any calibration, the proper gas must be flowing through the analyzer, some runs must be done and the readings must be stabilized.

To perform a calibration, it is suggested to perform the following steps:

- Configure your system (peak data, valve timings, system configuration data, alarm data, range data, flow tables and PID settings) and your calibration data (zero and span gas).
- Enable Noise Threshold Calibration with the corresponding button in the **CALIBRATION-CALIBRATION MENU**.
- Start a calibration with the start button. Ranges #1 will be set automatically.
- After some cycles have been done (you are supposed to see a relatively straight line in the RUN-REAL TIME CHROMATOGRAM MENU), press Calculate Threshold.
- Disable Noise Threshold Calibration with the corresponding button.
- Enable Span calibration, select impurity values and press Start.
- After some cycles have been done (you are supposed to see peaks), press **ReSpan**.
- Disable span calibration with the corresponding button.
- Your system is calibrated.



Calculate Threshold, ReSpan and Start buttons are only accessible when calibration is enabled.

The other values displayed on the grid are:

- Area counts: this is the area that covers the peak in counts (digital converter units). This area is used to generate the concentration displayed in the Result column;
- Response factor: this is the area counts divided by the value in the Result column.

Moreover, there is an **Automatic Re-Span** option. When this option is checked, the analyzer automatically performs a re-span (same as pressing the **Re-Span** button) after a cycle of calibration reaches an end. In other words, after having press the **Start** button in the **Span** calibration, the cycle starts, then ends and the analyzer automatically calculates the value of this calibration with the last measured area counts and displays it in the **Result** column.

#### 6.2.4.2 Calibration of the 4-20mA outputs

By clicking **Calibration** on the menu bar and then **4-20mA outputs**, you access the **4-20MA OUTPUT CALIBRATION MENU** (fig. 37).

- 4-20 ma Calibration
IO Board #1
4-20 mA #1
4-20 ma Output % (0.000 to 100.000): 78.996 Set
Set as low value MA
Set as high value mA
Calibrate

Figure 41 : 4-20 mA outputs calibration menu

This menu is used to calibrate the 4-20mA outputs. This is originally performed in factory. Therefore, you normally don't have to do it again.

- Step 1: First choose the IO board on which you want to calibrate the 4-20mA outputs; normally, the 4-20mA outputs are on the first IO board;
- ► Step 2: Select the output you want to calibrate;
- ► Step 3: Connect an ampere-meter on the rear panel of the analyzer on the corresponding output;
- ► Step 4: Output a value between 0 and 100% (advise: 20%) by writing it in the edit box on the left of the "Set" button and click on "Set";
- ► Step 5: Write in the edit box on the right of "Set as low value" the reading on your ampere-meter and press "Set as low value";
- ► Step 6: Redo step 4 and 5 with a value higher than the first (advise: 80%) but enter it in the "Set as high value" edit box;
- **Step 7**: Press on the Calibrate button.

### 6.2.5 Remote

#### 6.2.5.1 Report Software

Clicking on **Remote** on the menu bar and then on **Report Software** brings you to the **REMOTE-REPORT SOFTWARE MENU** (fig. 38).

	Address type • DHCP O Static
	IP Address : \$\$\$\$\$\$\$\$\$\$\$\$\$\$\$\$\$\$\$\$\$\$\$\$\$\$\$\$\$\$\$\$\$\$
	Subash Made 255 - 255 - 255 - 0
	Subnet Mask : 233 - 233 - 233 - 0
	Gateway : 192 . 168 . 75 . 1
	Set Static Parameters (E3)
	4
	Disconnect from Data Monitoring Software (F2)
	Test USB key for Report Software (F4)
4	

Figure 42 : Remote software menu

In this menu you can visualize and configure the IP address of the Servopro Chroma that you'll need to use in the Servopro Chroma Report Software.

Two type of IP address can be configured: Static or DHCP. A static address is an address that is manually set one time by the user of this analyzer. A DHCP address is an address that is automatically set by a router on the network. For more details, ask your Information Technology (IT) technician.

If you want to set a Static IP address, select the Static type, enter the three configuration parameters, then press the **Set Static Parameter (F3)** button.

The **Disconnect** button is there if you want to close the communication between the Servopro Chroma and the PC Software. <u>If you close the Servopro Chroma</u> <u>Report Software in the PC, you must use this button to tell the Servopro</u> <u>Chroma that the connection is terminated.</u>

Finally, the **Test USB key for Report Software** button is used to test if the analyzer detects a USB key that can then be used for acquiring data for the Report Software.

#### 6.2.5.2 Remote Control

Clicking on **Remote** on the menu bar and then on **Remote Control** brings you to the **REMOTE-REMOTE CONTROL MENU.** 

Remote Control Settings     Port :   1076     Password :   10006

Figure 43 : Remote control menu

In this menu, you are able to set the network port number on which the remote control software will communicate with the Servopro Chroma and the password used for the login connection.

### 6.2.6 About

Clicking on About on the menu bar brings you to the ABOUT MENU.

bout K4000	
K4000 Softwa	are Version 2.5.6.3
Build Time : 0	08/01/07 23:01:19
Contrôle Analytic	ue Copyright (C) 2005
www.cai-ca.com for more info	rmation on Contrôle Analytique Inc.
Module name	Firmware version
Conditionning board #1	1.52
IO Board board #1	1.34 / 1.14
Pressure board #1	1.32
Pressure board #2	1.32
Pressure board #3	1.32
Memory	
	-
Free disk space : 220 Mi	5
Free disk space : 220 Mi Free ram space : 69 MB	ок

Figure 44 : About menu

In this menu, you have the software version of the user interface (PC), the firmware versions of the embedded electronic boards and information about the memory usage. This information can be very useful if you have to contact Servomex support.

### 6.2.7 Chromatogram saving

In version 2.6.3, a new feature was added to store chromatogram on the internal hard drive. This feature can be configured from menu Historic/Chromatogram. To enable this feature, the check box must be selected. When selected, chromatogram will be saved at the end of each cycle. If not selected, chromatogram will not be saved.

K4000 - Chromatograms Historic								
Run	Diagnostic	Configuration	Historic	Calibration	Remote	About	Quit	
			I	✓ Enable Cl	hromatogra	ams Stori	ng	
		Percer	itage of al	lowed space	for chroma	atograms	storage used :	
								0%
		Delete C	hromatogr	ams		Tran	isfer to USB stick	

Figure 45 : Chromatogram saving menu

A space of 100 Mb is configured on the hard drive. This allow around one month of continuous data. To visualize data, insert a USB key into the port in front of analyzer and press "Transfer to USB stick" button. If USB key is not detected, an error message will be displayed. Chromatograms can be deleted by pressing delete button. This will delete all chromatograms on the hard drive.

Once chromatograms are transferred to the USB stick, copy them on a PC computer and use the import feature of ReportSoftware to import them.

The progress bar indicates the actual used space on the hard drive.

### 6.2.8 Quit

It is important to always use the Quit button of the menu bar when the analyzer **must be turned off.** If you close the analyzer without quitting, some data could be corrupted. After a quit is done, wait for the black screen and then turn the analyzer power supply off with the switch on the back of the Servopro Chroma.

When clicking on the **Quit** button, the analyzer will ask you for the general password of this analyzer. This password is, by default, the serial number of the analyzer, but can be changed into the **CONFIGURATION-SYSTEM CONFIGURATION MENU.** 

# 7 INSTALLATION AND SETUP

## 7.1 System power up

If your system has slave chassis, they must be powered on before the master chassis. If your system has an external PC, start all chassis and finally power on the PC.

To power off the system, properly exit the PC software and you can then close chassis in any order.

## 7.2 Electrical installation

## 7.2.1 Electrical safety



#### WARNING

Ensure that the electrical installation of the analyser conforms to all applicable local and national electrical safety requirements.



### WARNING

Obey the safety instructions given below when you install the analyser; if you do not, the analyser warranty may be invalidated, the analyser may not operate correctly, or it may be damaged.



#### WARNING

All signal terminals are separated from the analyser mains circuits by reinforced insulation. The terminals must only be connected to circuits that are themselves separated from mains voltages by an isolation method that provides at least this level of protection.

#### The following warnings must also be considered:

- The electrical supply coupler or plug must be easily accessible for disconnection from the electrical supply.
- The electrical supply circuit must incorporate a suitable fuse or over-current protection device, set to or rated at no more than 10 A for each individual chassis of the system.
- Ensure that your electrical supply can provide the necessary maximum power consumption: refer to Section 2.1.
- All signal and electrical supply cables must be rated of temperature of 70 °C or higher.
- Ensure that the cables that you connect to the analyser are routed so that they do not present a trip hazard.

- When you carry out insulation testing, disconnect all cables from the analyser.
- Ensure the analyser is provided with a sound earth connection via the electrical supply plug.

### 7.2.2 Master chassis connection

The following drawing explains how to connect the different parts of a Servopro Chroma system that only have one slave chassis. The master chassis must be connected to a power source using a properly rated cable (see section 3.13). If an auxiliary oven is required for your application, connect the cable coming out of the auxiliary oven to the connector named "Auxiliary oven" (see section 1.3.1).





### 7.2.3 Slave chassis connection

The following drawing explains how to connect the different parts of a Servopro Chroma system that have master and slave chassis. The master and slave chassis must be connected to a power source using properly rated cables (see section 3.13). If an auxiliary oven is required for your application, connect the cable coming out of the auxiliary oven to the connector named "Auxiliary oven". In case you have two auxiliary ovens, make sure that you connect the right oven to the right chassis.

In order to operate, the master chassis must be able to communicate with the slave chassis. Connect both RS-485 ports using the high density DB-9 cable provided with the system (see section 1.3.1).



Figure 47 – Master and slave chassis electrical connections



WARNING

The units shall only be connected using the cables supplied by Servomex.

## 7.2.4 Stand alone PC with slave chassis

The Servopro Chroma can be supplied with a stand alone PC depending on the application. The stand alone PC must always be connected to the same protective earth as the slave chassis of the system

The following drawing explains how to connect the different parts of a Servopro Chroma system that have a stand alone PC and slave chassis. The stand alone PC and slave chassis must be connected to a power source using properly rated cables (see section 3.13). If an auxiliary oven is required for your application, connect the cable coming out of the auxiliary oven to the connector names "Auxiliary oven". In case you have two auxiliary ovens, make sure that you connect the right oven to the right chassis. In order to operate, the stand alone PC must be able to communicate with the slave chassis. Connect the port "To analyser" on the PC to the RS-232 port on slave unit 1. If you have a second slave in your system, connect port "Aux" on the PC to the RS-232 port on slave unit 2.

Both connections should use the high density DB-9 cable provided with the system (see section 1.3).



Figure 48 – External PC connections



WARNING

The units shall only be connected using the cables supplied by Servomex.

## 7.3 Connecting Servopro Chroma rear connector

## 7.3.1 Digital input

Servopro Chroma has one digital input that is used to start an analysis remotely. Connection between Servopro Chroma and Remote system is made with a shielded twisted cable. In case of fuse failure, change fuse F12, referring to section 7.3.1.

The shield of the digital input contacts cable must be terminated to the analyser on one the "E" ports of the rear connector.



Figure 49– Digital input connection



## 7.3.2 Digital output contacts

Servopro Chroma can have up to 12 digital outputs. All those outputs are dry contacts. No power is provided by those outputs. They all connected to a common port and fused. Connection between Servopro Chroma is made using a shielded multi core cable. In the case of fuse failure, change fuse F14, referring to section 7.3.1.

Port Name	Purpose
$\mathbf{E}$	Earth point for cable screening
R1 – R8	Range ID contacts
A1 – A2	Alarm indicators
ST	Status indicator
EX1	Calibration indicator or Cycle indicator

Table 24 : Digital Output Contacts

The shield of the digital output contacts cable must be terminated to the analyser on one the "E" ports of the rear connector.



Figure 50 – Digital outputs connection



#### WARNING

The digital output terminals are separated from the analyser mains circuits by reinforced insulation. The terminals must only be connected to circuits that are themselves separated from mains voltages by an isolation method that provides at least this level of protection.

## 7.3.3 4-20 mA outputs

Figure 44 represents the circuit of Servopro Chroma 4-20 mA. The connection between the Servopro Chroma and the monitoring system must be done using a shielded twisted cable. Connect the load resistor on the monitoring system. A differential input must be used on the data acquisition system to properly monitor the data. In case of fuse failure, refer to Table 1 and section 7.3.1.

The twisted pair must be terminated at a matched pair (+ -) of numbered output ports and the shield must be terminated to the analyser on one the "E" ports of the rear connector. Each numbered output is set to a particular impurity represented on the chromatogram and can be changed via the configuration menu.

The HR output is the high-resolution chromatogram, it is the current reading of the GC cell as it progresses through a cycle.



Table 25: 4-20 mA connections and associated fuses



#### WARNING

The analogue output terminals are separated from the analyser mains circuits by reinforced insulation. The terminals must only be connected to circuits that are themselves separated from mains voltages by an isolation method that provides at least this level of protection.

## 7.3.4 FID fuel shut-off relay/Hydrogen relay output



#### WARNING

A fuel shut-off valve must be connected to the Servopro Chroma if a FID detector is installed to prevent the accumulation of fuel. The inside control valve is not considered tight shut-off.

A fuel shut-off valve can be connected to the rear connector of the analyser. This valve must be normally closed. This valve is controlled by the Servopro Chroma software. When an error occurs, it closes this valve automatically to prevent the accumulation of fuel inside the Servopro Chroma.

See the FID Addendum (Section 10) for more information when using a FID detector.



Figure 52 – Connecting a fuel shut-off valve



#### WARNING

The FID relay output terminals are separated from the analyser mains circuits by reinforced insulation. The terminals must only be connected to circuits that are themselves separated from mains voltages by an isolation method that provides at least this level of protection.

## 7.3.5 RS-485 port

The Servopro Chroma has a RS-485 port. This port is only used to communicate with other Servopro Chroma systems. It is not design to be used to interface with third party systems. If your system has multiple chassis, you need to connect them using shielded D-sub HD15 Male/Male cable.



#### WARNING

The RS-485 output terminals are separated from the analyser mains circuits by reinforced insulation. The terminals must only be connected to circuits that are themselves separated from mains voltages by an isolation method that provides at least this level of protection.

### CAUTION

To comply with EMC requirements, the RS-485 shall only be connected using the supplied cable. Refer to section 7.3.2 before making any electrical connections.

## 7.3.6 RS-232

With this option installed, this analyser can be remotely controlled using a RS-232 port. The computer is connected to the analyser by the mean of a DB-9 RS-232C (Null-modem) cable. The computer must have appropriate software to read the data transmitted by the analyser. See the RS-232 addendum for information about the analyser communication protocol.



### WARNING

The RS-232 output terminals are separated from the analyser mains circuits by reinforced insulation. The terminals must only be connected to circuits that are themselves separated from mains voltages by an isolation method that provides at least this level of protection.

## 7.3.7 Auxiliary oven port

The Servopro Chroma system can use auxiliary oven to perform some applications. The auxiliary oven normally contains a trap. The auxiliary oven must be connected to the auxiliary oven port on the rear panel. No third party oven can be connected to this port.

## 7.4 Gas installation



### WARNING

It is of the highest importance to never pressurize the analyzer because the quartz cell would be irreversibly damaged. So, before supplying any gas to the instrument, first <u>remove</u> the caps on the <u>vent</u> connections. However, in such a way as to not contaminate the analyzer it is better to leave all caps on gas connections until the gas installation is done and properly purged.



#### WARNING

This analyser is not to be used with any sample gasses that form a flammable mixture.

## 7.4.1 Introduction

The sampling system is the most important part of your analytical system. The performance of your analyzer can be dramatically limited by your gas transporting system. What we mean, by gas transporting system, any pressure regulator, valve, line, fitting, filter, purifier, etc. which are in contact with the sample or carrier gas to be introduced into the analyzer.

Absolutely, no leaks can be tolerated. Here, by a leak, we mean introduction of outside contaminants originally not present in the system. In our case, such leaks are atmospheric inboard leakage.

For all gas line connections (including calibration and sample), we recommend the use of 1/8" stainless steel tubing in full length, no fittings. You must avoid pipe thread connections, because they are usually sealed with Teflon tape and some particles can be introduced into the lines. By using compression type tube fittings, the venturi aspiration of outside contaminants is virtually eliminated.

Tube quality is often overlooked. For 1/8" O.D. tube, use a minimum wall thickness of .028". The tube must be purchased to meet ASTM69 specifications. Inferior quality tubes may have irregularities on their circumference causing bad sealing with compression tube fittings.

Install a by-pass rotometer on your sample line. It must be installed near the analyzer and close to the stream selection valve. This increases the gas velocity and allows a faster purging of the sample line, before selection. You will also get a faster response time this way. Also adjust your sample line pressure to a value, which will bring proper flow in the system. All lines must be cleaned and purged to remove any traces of moisture or particles. Particles can damage your stream selection valves. Also, particles in the inlet filter will trap moisture.



The best sampling system available today is our MSS, i.e. Manual Sampling System or our ISS. These will provide many years of trouble free service. They are easily remotely controlled, perfect for Auto Calibration Systems, etc. Please contact us or visit our Website for more information.

### 7.4.2 Carrier Gas Hardware Selection

We strongly recommend the use of a gas purifier fed by UHP grade gas cylinder as a source of carrier gas. The purer the carrier gas, the better the performance of the analyzer. The cost associated with research grade cylinder of argon is high. The carrier grade argon cylinder quality is not enough.

The least costly solution over time, and better for the quality of carrier gas is to use a heated type getter based purifier. The cost of UHP grade argon cylinders is inexpensive compared to carrier or research grade. The purifier will last many years.

#### Gas Purifier

You can buy such purifier from Servomex's representatives

The model number to order is:	GP-200- 120: 120 VAC
	GP-200- 240: 240 VAC

You must add the proper model definition digits for supply voltage and fittings type. **See purifier specification sheet.** 

#### Cylinder Pressure Regulator

The pressure regulator must be a double stage type and made of stainless steel. The outlet pressure range must be selected according to the carrier pressure setting indicated in the configuration in the back cover of this manual.

The pressure stability is a critical parameter affecting the accuracy of the analyzer. A good quality pressure regulator must be used. For calibration gas cylinder, an outlet pressure of 200 kPag (30 psig) maximum is recommended.

#### Isolation and Samples Stream Selection Valves

The valves used for isolation or sample selection must be packless type i.e. diaphragm or bellows type. This is the only way to remove air diffusion.

## 7.4.3 Connecting the Servopro Chroma

- 1. Once Servopro Chroma installed in a cabinet or rack, remove all caps. It is very important because avoiding removing caps on the outlet will damage the plasma cell.
- 2. Connect all carrier gas inlets. Make sure that pressure are according to specifications and that lines were properly purged. Up to two carrier gasses may be used.
- 3. If your system has a trap or an auxiliary oven, install it. Connect the trap or oven inlet first.
- 4. Connect sample lines. Make sure that pressures are set as specified in the specifications. Up to two sample gasses may be used.
- 5. Connect all vents. These lines should be unpressurised, as pressurising the detectors within the Servopro Chroma will damage them.
- 6. Power on the analyzer and wait for stabilization.

# 8 GENERAL OPERATION

### 8.1 Routine maintenance

The analyzer does not need a lot of supervision. But for accurate measurements you must be sure that:

- 1. Carrier flow is as per the value specified for your configuration. To check the carrier flow, put the injection mode in manual. Go back in «RUN» menu and watch the value. It is normal to have carrier flow deviation during a cycle. Between and before an injection the carrier flow must be correct.
- 2. From time to time, verify the oven temperature in the diagnostic menu. It must be constant and at the value specified in the configuration section in the back cover.
- 3. The sample flow must be at the same value of the previous calibration procedure. This is to make sure you have the same sample volume injected. The default value of 75 cc/min is OK.
- 4. Clean cabinet fan's filter periodically. This is very important, high internal temperature will eventually cause problems.

# 9 STORAGE AND DISPOSAL

## 9.1 Storage

Refit any protective cover (see Section 5) and place the Servopro Chroma and any associated equipment in its original packing before storage.

Store the Servopro Chroma and any associated equipment in a clean, dry area. Do not subject it to excessively hot, cold, or humid conditions: See section 2.5.

# 9.2 Disposal

Dispose of the Servopro Chroma and any associated equipment safely, and in accordance with all of your local and national safety and environmental requirements.



The Servopro Chroma is not suitable for disposal in municipal waste streams (such as landfill sites, domestic recycling centres and so on). Refer to section 13 for disposal requirements in accordance with the WEEE Directive within the EC.

# **10 COMPLIANCE**

- The analyzer complies with the European Community "Electromagnetic Compatibility Directive":
  - Emissions: Class B Equipment suitable for use in domestic establishments and in establishments directly connected to a low voltage supply which supplies buildings for domestic purposes.
  - Immunity: "Basic" Considered appropriate to equipment intended for use in domestic, commercial and light industrial environments.
- The analyzer complies with the European Community "Low Voltage Directive", by the application of EN 61010-1 and rated for Over Voltage Category II, Pollution Degree 2.
- The analyzer complies with the Class B digital apparatus requirements of ICES-003 of Canada through the application of EN 61000-6-3:2007.
- L'analyseur est conforme aux Conditions B numériques d'appareillage de classe de NMB-003 du Canada par l'application du EN 61000-6-3:2007.
- This analyzer complies with Part 15 of the US FCC Rules for Class B equipment. It is suitable for operation when connected to a public utility power supply that also supplies residential environments.
- The analyzer has been assessed to IEC 61010-1 for electrical safety including any additional requirements for US and Canadian national differences.

# **11 FID ADDENDUM**

## 11.1 Introduction

This document explains how to set up and use a FID with the Servopro Chroma software.

## 11.2 Overview

The low noise flame ionization detector (FID) is designed to measure hydrocarbons in the sub ppm level to percentage level. Measurement of atmospheric air for low level hydrocarbons and hydrocarbons content of exhaust emission from internal combustion engines are typical applications.

## 11.3 Logarithm based amplifier

Contrôle Analytique has designed a circuit allowing the measurement of weak currents in order to interface some of its detectors. Other analyzers based on FID detector use electrometers based on high resistance values in order to amplify the current. However, it is difficult to obtain good resistances in this type of circuit, which produces variations with fluctuations in temperature. Moreover, these resistances introduce noise, according to Johnson's Law, which states that noise caused by a resistance is proportional to its value and temperature. Contrôle Analytique therefore became innovative by making use of the perfect logarithm feature of transistors to design a logarithm based electrometer that is not only extremely stable with variations in temperature but also very quiet. The use of transistors has thus allowed us to eliminate the high value resistances, but also the noise and drift that they cause. Also, this circuit has made it possible to measure a vast range of currents, from femtoampere (10⁻¹⁵) to the milliampere (10⁻³). It has greatly improved the quality of the measurement made with Contrôle Analytique's detectors.

## 11.4 **MENUS**

### 11.4.1 CONFIGURATION-ADVANCE MENU



You normally shouldn't have to change these configurations; they are set in factory and reflect the hardware of the analyzer.

In the menu Advanced Configuration, you can add or remove FIDs by modifying the quantity.

10	1 <mark>0 - Advanc</mark> Diagnostic	ed Configu	uration tion Historic	Calibration	Remote	About	Quit				
	Diagnosae	Coningerer	don inscore	Calibration	Kelhote	MDOOL	Zaic				
G	eneral										
D	etector			General							
	1										
				Item	Qu	antity	- System Alarm				
FI	low			Cond. E	oard 1		⊙ On C Off				
Va	alve			FID I/O Boa	1 rd 1	2	RS-232 Communication				
0	ven			Flow	1		O On 💿 Off				
F	iD			Valve	1						
				Detecto	nr 2		Take Snapshot				
				Oven	3						
				Plasma	1						
				Press. B	Board 1						
				Temp. (	Itrl. O						
				D/O Boa	ard O						

Figure 53: Advanced menu

## 11.4.2 CONFIGURATION-ADVANCE FID MENU

In this menu you will be able to modify the FID settings. You have to click on save for your modifications to take effect.

K4000 - Advanced Configuration										
Run	Diagnostic	Configur	ation Histo	ric Calibrat	ion Remote	About (	Quit			
G	eneral etector 1 2					F	ID			
			FID #	Polariz.	Pressure	Carrier*	Air*	Fuel*	Steps	Seconds
E	low		FID #1	93.00	392.40	50.00	500.00	30.00	5	10
0	ven ID				*	Those set-p	oints will be t	used when th	ne flame mus	t be started
				ļ	Sav	e		ancel		

Figure 54: Advanced FID menu

Parameter	Range	Unit	Description					
Polariz.	0-200	Volt	Voltage that is applied to the FID polarizer.					
Pressure	300-500	In of H ₂ O	This value is used to calibrate the pressure sensor used by the FID. When save is pressed, the pressure sensor is calibrated with the value entered in the box.					
Carrier*	0-100	cc/min	This carrier set point is only used to ignite the FID. This set point is not used during normal operation.					
Air*	0-1000	cc/min	This air set point is only used to ignite the FID. This set point is not used during normal operation.					
Fuel*	0-200	cc/min	This fuel set point is only used to ignite the FID. This set point is not used during normal operation.					
Steps	0-10	Steps	Number of steps for ramping flow to normal value					
Seconds	0-60	Seconds	Time between each step.					
Table 26 : Advanced FID menu parameters								


Air, Fuel and carrier set points are for the ignition, if you press ignite flame in the Diagnostic menu, the Servopro Chroma will automatically change these set points to ignite flame and it will set back the normal value when the flame is lit.



## 11.4.3 CONFIGURATION-ADVANCE DETECTOR MENU

n C	Diagnostic	Configuration	Historic	Calibration	Remote	About	Quit			
Ger Det	ector				[	)ete	ctor #′	1		
Flov Valv Ove FID	w ve :n				FID	tor Type		•		
		C 1	iond. Board	i# (	Channel # I		Starting Co	punt	Location FID #1	<b>_</b>
					Save			ancel		

Figure 55: Advanced detector menu

In this menu you will be able to modify the FID detector settings. You have to click on save for your modifications to take effect.

- Cond. Board: Select on which acquisition board the FID is connected.
- Channel #: Select on which channel the Fid is connected on the acquisition board.
- Starting Count: Threshold value read by the detector inside the FID to determine if the flame is lit.
- Location: The position of the detector.



## 11.4.4 CONFIGURATION-ADVANCE FLOW MENU

This menu gives you the possibility to give a name to each flow associated with the FID. To operate, a FID needs Air, Fuel and Carrier gas. So, pressure regulators associate to these gases must be configured. A name must be given to each one. To set a name for Air, select Air in the location field and then enter the name you want into the field Name. Make sure to select the right Parent and the FID type. You must do the same for Fuel and Carrier.

General	_							
				Flow				
Jetector								
-								
Z								
low	Name	Parent	_	Location	_	Туре	Plasma	
alve	Sample	Regulator #1	•	IO Board #1	-	Sample 💌		-
	Carrier	FID #1	•	Carrier	•	FID 🔽		-
Jven	Air	FID #1	•	Air	•	FID 🔽	1	•
D	Fuel	FID #1	•	Fuel	•	FID 💽		-
		Sav	e		_			

Figure 56: Advanced flow menu



## 11.4.5 CONFIGURATION-ADVANCE OVEN MENU

The oven menu also works the same way. You can now select FID as a location, and you will be forced to use Channel #1 as the FID only allows configuring its first oven.

K40	00 - Advance	ed Configurati	on						
Rur	n Diagnostic	Configuration	Historic Calibral	tion Remote	About	Quit			
	General Detector				о	ven			
	1 2								
F	low		Name	L	ocation		Channel #		
	(ali ia		Oven 1	I/	O Board	#1	Channel #1	-	
			Oven 2	I/	O Board	#1	Channel #2	-	
	Oven		Oven FID	F	ID #1		Channel #1	-	
F	FID			Save			Cancel		

Figure 57: Advanced oven menu



## 11.4.6 MONITORING AND CONFIGURATION MENUS

Throughout the application, the flows and ovens associated with your FID will be showed with the other flows and ovens in your analyzer and you will be able to interact with them in the same way you would interact with the others. The 5 next figures will show you some examples.

Inside the System Configuration menu:

K4000 - System Configuration			
Run Diagnostic Configuration H	Historic Calibration Remote	About Quit	
Analog output Analog output Off C Track C Hold C Low C High	e Mode Generator Sta	rting Mode Range Mode Automatic	Auto-Start on Boot-Up C Yes © No
Flow Set Point Flow Name CC/min	Value CC/min	Oven Set Point Set Point Oven °C	Value °C
(Fuel) 0.0	0.0	Oven 1         50.0           Oven 2         120.0           Oven FID         0.0	1.8 0.4 29.4
Remote Starting Enable	G00 Sec	Range Configuration Active State Peak # on Range #2	Lock Range (F2)
Reload Manufacturing Setting	Print configuration	1 Opened 2 Opened	Set all opened
Change Password	Active Config : None		Set all closed

Figure 58: Monitoring fig 1

Inside the Real Time Chromatogram menu:

<b>K</b> 4	000	- Real T	ime Chro	matog	gram							
Ru	JN	Diagnosti	.c Configu	ration	Historic	Calibration	Remote	About	Quit			
	Al	Name	Value	R	All Acq. B	oard Acq.	Plasma 1			10/	11/2007 / 6:0	09:39
		1	< 1.00	1	— Ac	:q. #1	<ul> <li>Proces</li> </ul>	s Peaks				
		2	99999.00	2								
	Flow	/s										
		Name	CC/min									
	Air		0									
	Fu	el	0									
	Sa	mple	0									
	$\mathbb{P}$	<u> </u>										
	<u>ر</u>	ven										
		Name	/ °C			4.					0 / 600	sec
	E	Oven 1	0.5		Start (F2)	Historic	Values (F3	() His	toric Alarms (F4)	Range (F5)	Undo Zoom	(F8)
	Ľ	Jven 2	28.6	E F	Injection S	itatus	Inje	ction Mod	ie			
	1	10	20.0		On	C Off	0	Automat	ic 💿 Manual	Save (F6)		-2)
										Remote Cou	ntdown :	1 sec

Figure 59: Monitoring fig 2

Inside the Trending menu:

K400	0 - Trendin	g									
Run	Diagnostic	Configuration	Historic	Calibration	Remote	About	Quit				
						7					
4											F
۶.	art (F2)	enerator Power	(F3) R	e-Zero (F4)	R	.e-Zero D	et. (F5)			Time : 0 / 199	
A	ctive Detecto	r 🛛 Manual Valve	e   D/O Bo	ards	-1	Print (	Alt-F2)	sa sa	ave (Alt-F3	i) Undo:	Zoom (F8)
	Det.#	Acq # Gair	Fact	or Preamp	Flow	s		- 1		Oven	
11		1 0		0	Nam	e (	IC S	5etPnt	Follow	Name	°C
ŀ	2	1 0	x0	0	Air		) (	)		Oven 1	0.3
					Fuel		) (	)		Oven 2	0.0
					Sam		, L	J			28.4
								_			

Figure 60: Monitoring fig 3

Inside the PID Configuration menu:

K400	0 - PID Con	figuration					
Run	Diagnostic	Configuration	Historic	Calibration	Remote	About	Quit
	Flo	W Air Proportional Integrator Derivative	2.00			-Ov	en FID Proportional 5.00 Integrator 0.02 Derivative 0.00
			iave				Save

Figure 61: Monitoring fig 4

0 in	<b>.</b>	Control pop disa	ole					
AI								
Flow	PSI							
0.0	0.0							
0.0	0.0							
0.0	0.0	Flow Settin	ig tools					
0.0	0.0				_			
0.0	0.0			0.0	0	PSI		
0.0	0.0						Flow C	Control Type
0.0	0.0						C	Direct
0.0	0.0			◀	•			Direct
0.0	0.0						C	Bypass
0.0	0.0					1		
0.0	0.0		PSI :	2.3				
0.0	0.0							
0.0	0.0							
0.0	0.0	Actual co	ntrol outpu	it:   100.	0%			
0.0	0.0	Flow	value (CC)	): 0.0				
0.0	0.0			,				
,		Eactor :		Slop	🗖			
	- (L)			Diop	·			

Inside the Flow Configuration menu:

Figure 62: Monitoring

## 11.4.7 DIAGNOSTIC-DIAGNOSTIC MENU

In the Diagnostic menu, you will be able to change the FID parameters.

You'll have to click on the FID Board #X tab and then you'll be able to set its Gain, Offset and Pre-Amp. <u>These values are used in the trending menu.</u> You can also use these values to configure the default values in the Configuration/Cycle menu. You will also see its internal temperature as well as the flame's status. Finally, you'll be able to manually ignite the flame by clicking on the *Ignite Flame* button.

IN	Diagnostic	Conf	iguration	Histor	ric Calibration	n Remote	About	Quit				
Ac	q. Board #1											
[	-Detector si	gnal —				Gain:	7		Scale F	actor: 1	×	-
			Count		Volt		<u> </u>				· ·	
	Cell		375836		0.112008	Pre-Amp:	1	-	Po	larity: F	ollower	•
	Chromato	gram	8389127		2.500155							
	Detector #			•	1				Zero (F8)			
10	Board #1	FID Bo amplij	oard #1					- Pressure F	Regulator			
	Gain :		0.00		Official 1	E 050000		Flow	Count	CC/min	PSI	%
	Gailt		0.00		AISEC.	5.050000		Sample	584	0.0	1.0	0.0
	Pre-Amp :	8	0					Carrier	18472	15.0	28.0	69.0
		,t —					- 11	Air	579	0.0	0.9	83.1
	-Environmen			Due	<i></i>	444.6		Fuel	612	0.0	0.9	42.1
	-Environmen Temp. (°C)	: [9	11.6	Pre	ss. (inH2O) :	1					1	

Figure 63 : Diagnostic menu

## 11.4.8 CONFIGURATION-CYCLE MENU

In the Cycle menu, you will be able to change the Log Amplifier parameters that are associated to each FID peak.

When you have right selected the parameters in the Diagnostic menu, you have to set them in the cycle menu to be sure these parameters will be set on each cycle. To do so, go into peak/Configuration and select a FID peak. The peak data panel will allow you to enter FID Boards parameters like gain, offset and Pre-Amplification.

General	Name	Avr	Start	End	Det	Alarm 1	Alarm 2	Range 2	Range 1	Unit
Peak	C4	1	160.0	212.4	1 🔻	1000000.0	1000000.0	5	1x 💌	PPM
Configuration	C4H6	1	220.0	230.0	1 💌	1000000.0	1000000.0	5	1x 💌	PPM
Detection	CH4	1	264.4	305.0	1 💌	1000000.0	1000000.0	100	1x 💌	PPM
	C2H6	1	329.6	391.2	1 💌	1000000.0	1000000.0	5	1× 💌	PPM
Valves	C2H4	1	410.0	500.0	1 💌	1000000.0	1000000.0	5	1x 💌	PPM
Flow	C3H8	1	598.0	754.0	1 💌	1000000.0	1000000.0	5	1x 💌	PPM
Scripts	C2H2	1	781 8	930 N	1 🔻	1000000.0	1000000.0	5	1v 🔻	
	-Peak (	¢1 da	ta							_
										02
	Co	nd. B	oard :	Gain:	13		Plasma	power:	0	70
	Co	nd. B	oard :	Gain: Polarity:	13 Follow	er 💌	Plasma Pre-Amp	power: lification:	0	•
	Cc F1	ind. B D Boa	oard : ard :	Gain: Polarity: Gain:	13 Followe	er 💌	Plasma Pre-Amp	power: lification: Offset:	0 1 5.0100	•

Figure 64: Cycle menu

#### 11.5 Startup

- 1. Connect all gases to the analyzer and make sure that the tubes were properly purged.
- 2. Turn on the analyzer
- 3. Set flow set points to the following values for purging:

Gas	Flow set point (cc/min)
Fuel	0
Air	300
Carrier	50
Table 27 : Gas flow set point	

4. Wait for FID temperature stabilisation. Once temperature is stable, set flow set points to the following values:

Gas	Flow set point (cc/min)
Fuel	40
Air	300
Carrier	50
Table 28 : Gas flow set point 2	

Note: Those data can be different. See document settings that were sent with your Servopro Chroma.

5. Once flows are stable, press the ignite button to start the detector. The flow set point will automatically change to ignition set point and the detector will ignite.

Gas	Flow set point (cc/min)
Fuel	80
Air	500
Carrier	50
Table 29 : Gas ignition Point	

Note: Those data can be different. See document settings that were sent with your Servopro Chroma.

- 6. When the flame is lit, the flow set points will come back to there normal values. (Step 4)
- 7. The Servopro Chroma /FID is ready to use.

## 11.6 System errors

This section describes in details all possible system errors.

Air flow deviation: This error occurs when the difference between air flow and air flow set point is higher than 5 cc/min. When this error is active, it is impossible to ignite the flame.

Low air flow: This error occurs when the air flow is below 10 cc/min. When this error is active, it is impossible to ignite the flame.

**Fuel flow deviation:** This error occurs when the difference between fuel flow and fuel flow set point is higher than 5 cc/min. When this error is active, it is impossible to ignite the flame.

Low fuel flow: This error occurs when the fuel flow is below 10 cc/min. When this error is active, it is impossible to ignite the flame.

**Sample flow deviation:** This error occurs when the difference between sample flow and sample flow set point is higher than 5 cc/min. When this error is active, it is impossible to ignite the flame.

**Ambient temperature deviation:** This error occurs when ambient temperature is out of range. Ambient temperature must be between 5°C and 50°C.

**F.I.D. temperature deviation:** This error occurs when the difference between actual F.I.D temperature and F.I.D. temperature set point is higher then 10°C. When this error is active, it is impossible to ignite the flame.

**Glow plug damaged:** A glow plug is used to ignite the detector. If the glow plug is damaged, this error message is displayed. When this error is active, it is impossible to ignite the flame. The glow plug must be changed.

Flame off: This error occurs when the FID's flame is off.

**Sample flow reg. problem**: A problem appears with sample pressure regulator, the system will have to be restarted. If after rebooting, you still have this error, contact Servomex.

**Air flow reg. problem**: A problem appears with air pressure regulator, the system will have to be restarted. If after rebooting, you still have this error, contact Servomex.

**Fuel flow reg. problem**: A problem appears with fuel pressure regulator, the system will have to be restarted. If after rebooting, you still have this error, contact Servomex.

Every time an error occurs, the status contact relay is closed. Moreover, if you have set the "mA failure mode" to low or high, the 4-20 mA analogue output will be set according to this setting.

Because the SERVOPRO FID uses explosive gas, the software has built-in functions to reduce the risk of explosion. If the flame is off for more than 5 minutes and that the set point of hydrogen is not 0 cc/min, it will be automatically set to 0 cc/min and the hydrogen shut off valve will be de-activated.



When one of the gas is not flowing in the analyser, it is normal that the flame goes off. It is also normal that the flame goes off when there is no sample gas into the analyser.

# 11.7 Troubleshooting

Symptom	Problem cause	Possible remedy	
	Fuel gas emerging from burner jet diluted with other gases because fuel system is insufficiently purged	Purge fuel gas regulator.	
Burner doesn't ignite	Air and fuel flow not properly set.	Adjust air and flow set points.	
	Glow plug damaged. Analyser not able to reach starting flow set points	Install a new glow plug. Verify all inlet pressures.	
	Air or fuel is contaminated.	Replace air or fuel cylinder. Add hydrocarbon traps. Purge tubing	
	There is a leak on sample line.	Leak test your sampling system.	
Analyser's reading is noisy	Flows are not stable	verify that all gas pressures are stable. Verify that flow readings are stable. Optimize PID loop constants.	
	Tubing used to connect gases to the SERVOPRO FID was not properly purged.	Clean tubing	
	There is a dead volume on one of the gas stream.	Verify plumbing diagram and remove any dead volume.	
	Exhaust is blocked by water	Make sure the exhaust tubing has a slope.	
Loss of sensitivity or no	Electrical elements of burner partially short circuited by combustion products.	Clean burner.	
	Polarisation spring is not in contact with FID jet.	Open FID detector and verify that the spring is in contact with the jet.	
False flame indication	Flame sensor may be damaged.	Change flame sensor.	

 Table 30 : Troubleshooting

# 12 RS-232 ADDENDUM

The communication parameters of the analyser are:

- Baud rate: 38400
- Parity bits: none
- Data bits: 8
- Stop bits: 1

The following tables describe the command set:

### 12.1 Start analysis

					K4000	
Request / Command				R	esponse	
ASCII Code					ASCII Code	
Character No	Character	No (hex)	Character No	Character	No (hex)	Meaning
			1	>	3Eh	
			2	А	41h	command ID
1	>	3Eh	3	1	31h	Analysis started
2	А	41h	4 (checksum)	q	72h	
3 (checksum)	А	41h	5	<	3Ch	
4	<	3Ch	1	>	3Eh	
			2	А	41h	command ID
			3	0	30h	Failure
			4 (checksum)	r	71h	
			5	<	3Ch	

Table 31 : RS-232 Start Analysis

### 12.2 Analyser status

						K4000		
Requ	est / Comm	and	Response					
		ASCII Code	ASCII Code					
Character No	Character	No (hex)	Character No	Character	No (hex)	Meaning		
			1	>	3Eh			
			2	В	42h	command ID		
			3	1	31h	1 = success		
			4	1 or 0	31h or 30h	Analyser ready: 1=ready; 0=not ready (still running)		
			5	1 or 0	31h or 30h	Run Modus: 1=measurement; 0=not measurement		
1	>	3Eh	6	1 or 0	31h or 30h	Run Modus: 1=calibration; 0=not calibration		
2	В	42h	7	1 or 0	31h or 30h	General Alarm: 1=OK; 0=Alarme active		
3 (checksum)	В	42h	8	1 or 0	31h or 30h	Flow Alarm Sample: 1=OK; 0=Alarm active		
4	<	3Ch	9	1 or 0	31h or 30h	Flow Alarm Carrier: 1=OK; 0=Alarm active		
			10	1 or 0	31h or 30h	Detector Alarm (Plasma off?): 1=OK; 0=Alarm active		
-			11	1 or 0	31h or 30h	Calibration Alarm: 1=last cal. OK; 0=last cal. invalid		
			12	1 or 0	31h or 30h	Measurement Alarm 1=last measurement OK 0=invalid		
			13 (chksm)	Х				
			14	<	3Ch			
			1	>	3Eh			
			2	В	42h	command ID		
			3	0	30h	0 = failure		
			4 (chksm)	r	72h			
			5	<	3Ch			

#### Table 32 : RS-232 Analyser Status

## 12.3 Send results

	K4000 Response				nmand	uest / Con	Re
	Meaning	ASCII Code No (hex)	Character	No	ASCII Code No (hex)	Character	No
	Command ID	3Eh 2	č	1			
154 (checksum))*	0 = Failure, 1 = Success (M failure, jump to Char no 154 (checks) Examples of format of Concentration Result from Pask No. 1: 13.00 (could be 15.00 %, 15.00 gpm or 15.00 ppb) 1000000.00 (could be 1 Mio. ppm or 1 Mio.ppb)	30h or 31h 20h or 30h to 39h 20h or 30h to 39h 2Eh	D or 1 Space or any Character between 0 and 9 Space or any Character between 0 and 9	3 5 7 8 9 10 11	3Ch 43h 43h 3Eh	> C C <	1 2 3 (chksm) 4
=ppm; 4=ppb	Semicolon for Data Separation Dimension of concentration of Peak No. 1: 1=%; 2=ppm; 4=pp Semicolon for Data Separation measuremerv subit for Peak No. 1: 1=valit; 0=invalid	20h or 30h to 39h 38h 31h or 32h or 34h 38h 31h or 30h	any Character between 0 and 9 : 1 or 2 or 4 : 1 or 0	13 14 15 16 17			
for Peak No. 2	Semicolon for Data Separation	3Bh 20h or 30h to 39h 20h or 30h to 39h	Space or any Character between 0 and 9 Space or any Character between 0 and 9	18 19 20 21 22 23 24 25 26			
=ppm; 4=ppb	Semicolon for Data Separation Dimension of concentration of Peak No. 2: 1=%; 2=ppm; 4=pp Semicolon for Data Separation	20h or 30h to 39h 20h or 30h to 39h 38h 31h or 32h or 34h 38h	any Character between 0 and 9 any Character between 0 and 9 : 1 or 2 or 4 : 4 or 0	27 28 29 30 31			
linvalid	measurement valid for Peak No. 2; 1+valid; 0-invalid Semicolon for Data Separation	31h or 30h 38h 20h or 30h to 39h 20h or 30h to 39h 20h or 30h to 39h	1 or 0 ; Space or any Character between 0 and 9 Space or any Character between 0 and 9 Space or any Character between 0 and 9	32 33 34 35 36			
tata for Peak No. 3	Concentration Result Peak No. 3:	20h or 30h to 39h 20h or 30h to 39h 20h or 30h to 39h 20h or 30h to 39h 20h or 30h to 39h 2Eh 20h or 30h to 39h	Space or any Character between 0 and 9 Space or any Character between 0 and 9 Space or any Character between 0 and 9 Space or any Character between 0 and 9 any Character between 0 and 9	37 38 39 40 41 42			
=ppm; 4=ppb	Semicolon for Data Separation Dimension of concentration of Peak No. 3: 1=%; 2=ppm; 4=pp Semicolon for Data Separation measurement valid for Peak No. 3; 1=valid; 0=invalid Semicolon for Data Separation	2un or 30h to 39h 38h 31h or 32h or 34h 38h 31h or 30h 38h	any unaracter between 0 and 9 : 1 or 2 or 4 : 1 or 0 :	43 44 45 46 47 48			
sta for Peak No. 4	Concentration Result Peak No. 4:	20h or 30h to 39h 20h or 30h to 39h	Space or any Character between 0 and 9 Space or any Character between 0 and 9 any Character between 0 and 9	49 50 51 52 53 54 55 56 57			
=ppm; 4=ppb	Semicolon for Data Separation Dimension of concentration of Peak No. 4: 1=%; 2=ppm; 4=pp Semicolon for Data Separation measurement valid for Peak No. 4: 1=valid; 0=invalid	20h or 30h to 39h 3Bh 31h or 32h or 34h 3Bh 31h or 30h	any Character between 0 and 9 : 1 or 2 or 4 : 1 or 0	58 59 60 61 62			
Data for Peak No. 5	Semicolm for Data Separation	38h 20h or 30h to 39h 20h or 30h to 39h 2Eh 20h or 30h to 39h 2Eh	Space or any Character between 0 and 9 Space or any Character between 0 and 9 any Character between 0 and 9 any Character between 0 and 9	63 64 65 66 67 68 69 70 71 71 72 73			
=ppm; 4=ppb	Semicolon for Data Separation Dimension of concentration of Peak No. 5: 1-1%; 2-ppm; 4-pp Semicolon for Data Separation measuremer valid for Peak No. 2; 1-valid; O-invalid Semicolon for Data Separation	3Bh 31h or 32h or 34h 3Bh 31h or 30h 3Bh 20h or 30h	: 1 or 2 or 4 : : 1 or 0 : : : :	74 75 76 77 78 79			
Data for Peak No. 6	Concentration Result Peak No. 6:	20h or 30h to 39h 20h or 30h to 39h	Space or any Character between 0 and 9 Space or any Character between 0 and 9 any Character between 0 and 9 any Character between 0 and 9	80 81 82 83 84 85 86 87 88			
=ppm; 4=ppb	Semicolon for Data Separation Dimension of concentration of Peak No. 6: 1–1%; 2–ppm; 4–pp Semicolon for Data Separation measurement valid for Peak No. 6; 1–valid; 0–invalid Semicolon for Data Separation	3Bh 31h or 32h or 34h 3Bh 31h or 30h 3Bh	1 or 2 or 4 1 or 0	89 90 91 92 93			
sk Døts for Pesk Nu. 7	Concentration Result Peak No. 7:	20h or 30h to 39h 20h or 30h to 39h 22h or 30h to 39h 20h or 30h to 39h 30h or 30h to 39h	Space or any Character between 0 and 9 Space or any Character between 0 and 9 any Character between 0 and 9 any Character between 0 and 9	94 95 96 97 98 99 100 101 102 103 104			
=ppm; 4=ppb a. ⊨invalid	Dimension of concentration of Peak No. 7: 1=%; 2=ppm; 4=pp Semicolon for Data Separation measurement valid for Peak No. 7; 1=valid; 0=invalid Semicolon for Data Separation	31h or 32h or 34h 3Bh 31h or 30h 3Bh	1 or 2 or 4 1 or 0	105 106 107 108			
Deak No. 8	Concentration Result Peak No. 8:	20h or 30h to 39h 20h or 30h to 39h 22h 20h or 30h to 39h 22h 20h or 30h to 39h	Space or any Character between 0 and 9 Space or any Character between 0 and 9 any Character between 0 and 9 any Character between 0 and 9	109 110 111 112 113 114 115 116 117 119			
-ppm; 4-ppb	Semicolon for Data Separation Dimension of concentration of Peak No. 8: 1=%; 2=ppm; 4=pp Semicoln for Data Separation measurement valid for Peak No. 8; 1=valid; O-invalid Semicoln for Data Secaration	38h 31h or 32h or 34h 38h 31h or 30h 38h	inty Character bereard of and 9	118 119 120 121 122 123			
or Peek No. 9	Concentration Result Peak No. 9:	20h or 30h to 39h 20h or 30h to 39h	Space or any Character between 0 and 9 Space or any Character between 0 and 9	124 125 126 127 128 129 130 131			
appm; 4=ppb d.	Semicolon for Data Separation Dimension of concentration of Peak No. 9: 1+W; 2+ppm; 4+pp Semicolon for Data Separation masarement valid for Peak No. 8; 1+valid; 0+irvalid Semicolon for Data Separation	20h or 30h to 39h 20h or 30h to 39h 38h 31h or 32h or 34h 38h 31h or 30h 38h	0 0 1 or 2 or 4 ; 1 or 0 ;	132 133 134 135 136 137 138			
its for Peak No. 10	Concentration Result Peak No. 10:	20h or 30h to 39h 20h or 30h to 39h 22h or 30h to 39h 22h or 30h to 39h	space or any Chanadre between 0 and 9 Space or any Chanadre between 0 and 9	139 140 141 142 143 144 145 146 147			
S-ppm; 4-ppb	Semicolon for Data Separation Dimension of concertration of Peak No. 10: 1-%; 2-ppm; 4-pp Semicolon for Data Separation measurement valid for Peak No. 10; 1-valid; 0-eivvalid Semicolon for Data Separation Checksum	2un dr sun to 39h 38h 31h or 32h or 34h 38h 31h or 30h 38h 32h	U : 1 or 2 or 4 : 1 or 0 : X ×	148 149 150 151 152 153 154 155			

 Table 33 : RS-232 Send Results

	Request / Command			Re	K4000 esponse	
Character		ASCII Code	Character		ASCII Code	
No	Character	No (hex)	No	Character	No (hex)	Meaning
1	>	3Eh	1	>	3Eh	
2	D	44h	2	D	44h	Command ID
3	0 or 1 (0 = disable $/ 1$ = enable)	30h or 31h	3	1	31h	Success
4 (chksm)	t or u	74h or 75h	4 (chksm)	u	75h	
5	<	3Ch	5	<	3Ch	
-			1	>	3Eh	
			2	D	44h	
			3	0	30h	Failure
			4 (chksm)	t	74h	
			5	<	3Ch	

#### Table 34 : RS-232 Enable-Disable Calibration

## 12.5 Respan

	Request / Command					K4000 Response
		ASCII Code			ASCII Code	
Character No	Character (meaning)	No (hex)	Character No	Character	No (hex)	Meaning
1	>	3Eh	1	>	3Eh	
2	E	45h	2	E	45h	command ID
3	Any number from 1 to 9 (number of the	31h to 39h	3	0 or 1	30h or 31h	0 = Failure, 1 = Success
	calibration point if multi-point)		4	u or v	75h or 76h	Checksum
4 (Peak 1)	0 or 1 (0 = don't respan, 1 = respan)	30h or 31h	5	<	3Ch	
5 (Peak 2)	0 or 1	30h or 31h				
6 (Peak 3)	0 or 1	30h or 31h				
7 (Peak 4)	0 or 1	30h or 31h				
8 (Peak 5)	0 or 1	30h or 31h				
9 (Peak 6)	0 or 1	30h or 31h				
10 (Peak 7)	0 or 1	30h or 31h				
11 (Peak 8)	0 or 1	30h or 31h				
12 (Peak 9)	0 or 1	30h or 31h				
13 (Peak 10)	0 or 1	30h or 31h				
14 (Chksm)	Х					
15	<	3Ch				

Table 35 : RS-232 Respan

## 13 Product disposal In accordance with the Waste Electrical and Electronic Equipment (WEEE) Directive

The label shown in Figure & is fitted to the analyser.



Figure 65 :The WEEE label

This label identifies that:

- The analyser is considered to be within the scope of the Waste Electrical and Electronic Equipment (WEEE) Directive.
- The analyser is not intended for disposal in a municipal waste stream, but shall be submitted for material recovery and recycling in accordance with the local regulations which implement the WEEE Directive.
- For additional information and advice on the disposal of the analyser in accordance with the requirements of the WEEE Directive, contact Servomex at info@servomex.com or your local Servomex agent.
- If you send the analyser to Servomex or your local Servomex agent for disposal, the analyser must be accompanied by a correctly completed decontamination certificate.

# **14 REACH Regulation**

In pursuance of the requirements included in Article 33 of the European REACH Regulation on the Registration, Evaluation, Authorisation and Restriction of Chemicals, information on Substances of Very High Concern (SVHC) contained in Servomex products is provided on <u>www.servomex.com</u>